X-RAY SOURCE

300 kV OPEN TYPE MICROFOCUS X-RAY SOURCE L12721

FEATURES

- ●High power: 300 kV
- No high voltage cable connection required
- Easy to replace cathode
- Easy operation
- High stability



SPECIFICATIONS

Parameter		Description / Value
Maximum tube voltage ^①		308 kV
X-ray tube voltage setting range		20 kV to 300 kV
X-ray tube current setting range	at 20 kV to 230 kV	0 μA to 1000 μA
	at 231 kV to 300 kV	0 μA to 500 μA
Cathode material		Tungsten
Minimum resolution		4 μm ^②
Maximum target power		Approx. 200 W
Target type		Reflection type
X-ray beam angle		40 degrees to 60 degrees
Target material		Tungsten
X-ray output window material / Thickness		Aluminum / 0.5 mm
Focus to object distance (FOD)		4.6 mm ± 0.4 mm
Protection		External short-circuit method (Normally closed)
Interlock 1 / Interlock 2		DC24 V / DC12 V
Input voltage (AC)	X-ray control unit	100 V to 240 V (50 Hz / 60 Hz) ^③
	Power supply for turbo pump controller	100 V to 240 V (50 Hz / 60 Hz) ^③
	Power supply for vacuum pump	100 V to 240 V (50 Hz / 60 Hz) ^③
Power consumption (Max.)	X-ray control unit	700 VA
	Turbo pump set	300 VA
Rated output		Continuous operation
Weight	X-ray tube unit / X-ray control unit	Approx. 111 kg / Approx. 11 kg
	Power supply for turbo pump controller	Approx. 1 kg
	Vacuum pump (diaphragm pump)	Approx. 4.3 kg
Operating ambient temperature	X-ray tube unit / X-ray control unit	+15 °C to +35 °C
Storage temperature	X-ray tube unit / X-ray control unit	+5 °C to +50 °C
Operating and storage humidity	X-ray tube unit / X-ray control unit	20 % to 85 % (No condensation)


NOTE: ①Maximum tube voltage during the warm-up period.

- 2X-ray chart use
- 3Auto switching




DIMENSIONAL OUTLINES (Unit: mm)

•X-RAY TUBE UNIT L12722

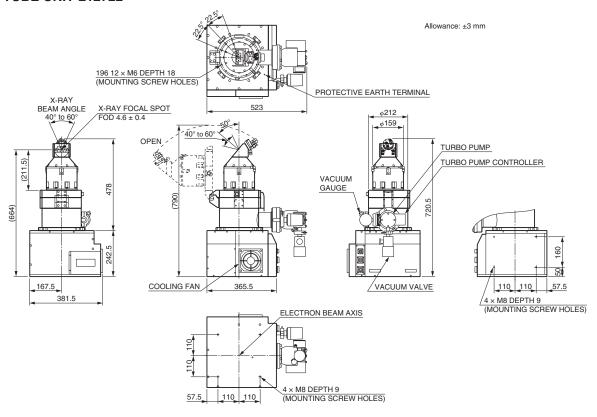

•X-RAY CONTROL UNIT C12723

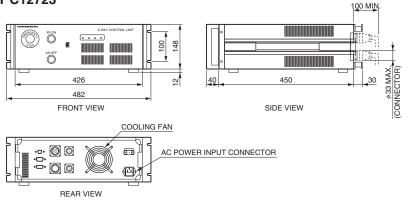

TXPRA0040EA

TXPRA0039EA

System configuration of L12721 X-ray tube unit (L12722) + X-ray control unit (C12723) + turbo pumps + accessories



PRECAUTIONS TO USE

· This microfocus X-ray source generates X-rays harmful to the human body. Use sufficient caution when handling the equipment to avoid direct or inadvertent exposure to X-rays.

Install the X-ray source or the X-ray tube unit in an X-ray shielded cabinet or room equipped with safety interlock functions to prevent accidental exposure to X-rays.

OPERATIONAL CAUTION

- · This microfocus X-ray source generates X-rays and must therefore be used only under the supervision of qualified personnel.
- This microfocus X-ray source shall be used in compliance with health and safety regulations enforced in order to prevent health hazards problems due to ionizing radiation.

Subject to local technical requirements and regulations, availability of products included in this promotional material may vary. Please consult with our sales office. Information furnished by HAMAMATSU is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications are subject to change without notice. No patent rights are granted to any of the circuits described herein. ©2019 Hamamatsu Photonics K.K.

HAMAMATSU PHOTONICS K.K. www.hamamatsu.com

HAMAMATSU PHOTONICS K.K., Electron Tube Division

314-5, Shimokanzo, Iwata City, Shizuoka Pref., 438-0193, Japan, Telephone: (81)539/62-5248, Fax: (81)539/62-2205

S14-5, Snimokanzo, Iwata City, Snizuoka Prer., 438-U193, Japan, Telephone: (81)539/62-5248, FaX: (81)539/62-22U5

U.S.A.: Hamamatsu Corporation: 360 Foothill Road, Bridgewater, NJ 08807, U.S.A., Telephone: (1)908-231-0960, Fax: (1)908-231-1218 E-mail: usa@hamamatsu.com

Germany: Hamamatsu Photonics Deutschland GmbH.: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49)8152-375-0, Fax: (49)8152-265-8 E-mail: info@hamamatsu.de

France: Hamamatsu Photonics France S.A.R.L.: 19, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 10 E-mail: info@hamamatsu.fr

United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10 Tewin Road, Welwyn Road, W

TXPR1032E01