

### ■ Features

- Installed MEMS grating
- Realize high speed and broad wavelength sweeping
- Built-in collimation lens
- Broadband QCL based on DAU structure


**Anticross  
DAU<sup>TM</sup>  
structure**



### ■ Outline

It is a pulse quantum cascade laser that realized broadly wavelength sweeping by external cavity configuration. It can carry out the measurement of mid-infrared spectrum remotely, contactless and with high throughput. This product offers new value to the mid-infrared spectroscopy, which was traditionally carried out with FT-IR, by utilizing features of laser; high output and high wavelength resolution.

### ■ External cavity

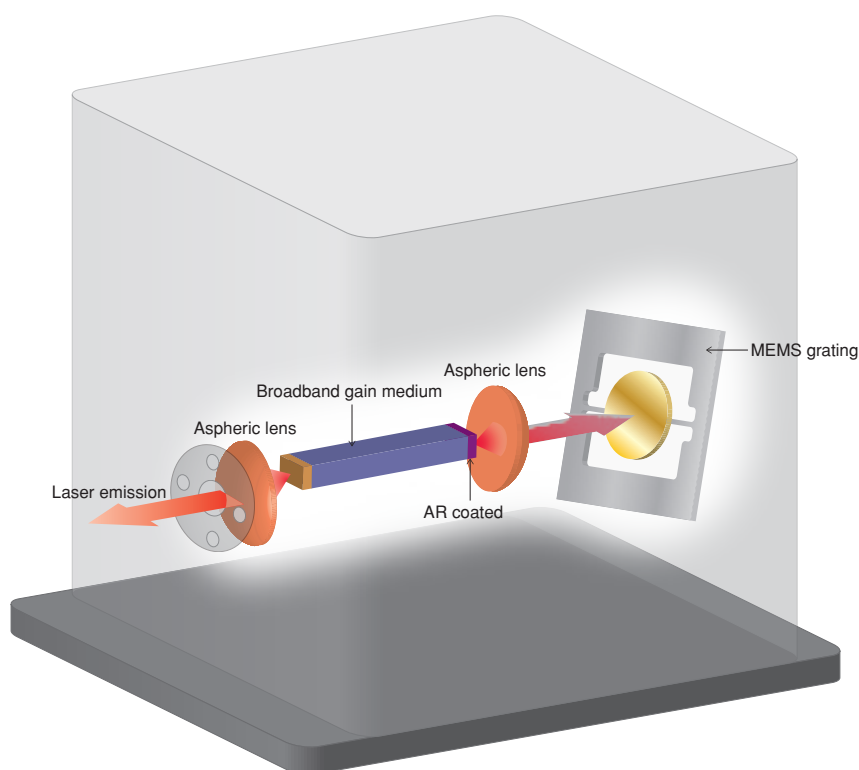

# Wavelength swept pulsed quantum cascade laser L14890-09

## ■ Characteristics

Unless otherwise specified, the repetition rate of the MEMS operation signals  $F(\text{mems})=1.8$  kHz, The repetition frequency  $F(\text{qcl})$  of the QCL operation signal is 180 kHz, and the QCL operation temperature  $T_{\text{op}(\text{qcl})}$  is 20 °C.

Characteristic	Condition	Min.	Typ.	Max.	Unit
Wavenumber sweep frequency	External control *1 (MEMS operation signal)	—	1.8 *2	—	kHz
Pulsed output power	MEMS stop	200	600	900	mW
Optical pulse width	Fixed before shipment from the factory	80	100	120	ns
Optical pulse repetition frequency	External control *1 (QCL operation signal)	—	180 *3	—	kHz
Spectrum linewidth *4	MEMS stop	—	2.0	3.0	cm <sup>-1</sup>
Beam diameter *5	Distance *6 800 mm ± 20 mm MEMS stop	3	5	7	mm
Beam divergency *7	Distance *6 800 mm ± 20 mm MEMS stop	—	—	5	mrاد
Beam steering stability	Distance *6 800 mm ± 20 mm	—	—	2	mrاد
Polarization	Vertical to the bottom				
Center wavenumber	Calculated from wavenumber sweep width	1020	1075	1140	cm <sup>-1</sup>
Wavenumber sweep width	Pulsed output power > 10 mW	180	200	—	cm <sup>-1</sup>

\*1: Controlled by a signal source such as a separate function generator.

\*2:  $F(\text{mems})$ : Repetition frequency of the MEMS operation signal.

\*3:  $F(\text{qcl})$ : Repetition frequency and rise edge of the QCL operation signal.

\*4: Full width at half maximum of the spectrum. \*5:  $1/e^2$

\*6: Distance from the outgoing window \*7: Half angle

## ■ Recommended operating conditions

Characteristic	Symbol	Value	Unit
Power input	Voltage	DC24	V
	Current	0.05	A
QCL operation signal *1	Waveform	Positive pulse square wave *2	—
	Repetition frequency	180	kHz
	Signalling voltage	4.0 *3	V <sub>pk</sub> *3
	Duty ratio	1.8	%
MEMS operation signals *4	Waveform	Sine wave	—
	Repetition frequency	1.8	kHz
	Signalling voltage	4.0	V <sub>pp</sub>
Peltier current *5	I <sub>c</sub>	±1.6	A
Peltier voltage *5	V <sub>c</sub>	±2.2	V
QCL operation temperature *5	T <sub>op(qcl)</sub>	+20	°C

\*1: External control. A signal source such as a separate function generator is needed. \*2: Positive pulse with a pulse base voltage of 0 V.

\*3: V<sub>pk</sub>: Pulse stop voltage value when the pulse base voltage is set to 0 V.

\*4: External control. A signal source such as a separate function generator that is different from \*1 is needed.

\*5: External control. Operation temperature of the laser element (QCL) controlled by a separate peltier controller.

## ■ General specifications


Characteristic	Value	Unit
Operation temperature (ambient)	+10 to +40	°C
Storage temperature (ambient)	+5 to +55	°C
Relative operation humidity (ambient)	30 to +60 *1	%
Relative storage humidity (ambient)	20 to +60 *1	%
Startup time	5 or less	Second
Place of use	Indoor place	—
Operation altitude	2000 or less	m
Dimensions (W × H × D) *2	82 × 88 × 112	mm
Weight	1.2	kg

\*1: No condensation \*2: Excluding projections

# Wavelength swept pulsed quantum cascade laser L14890-09

## ■ Configuration

### ● Connection example

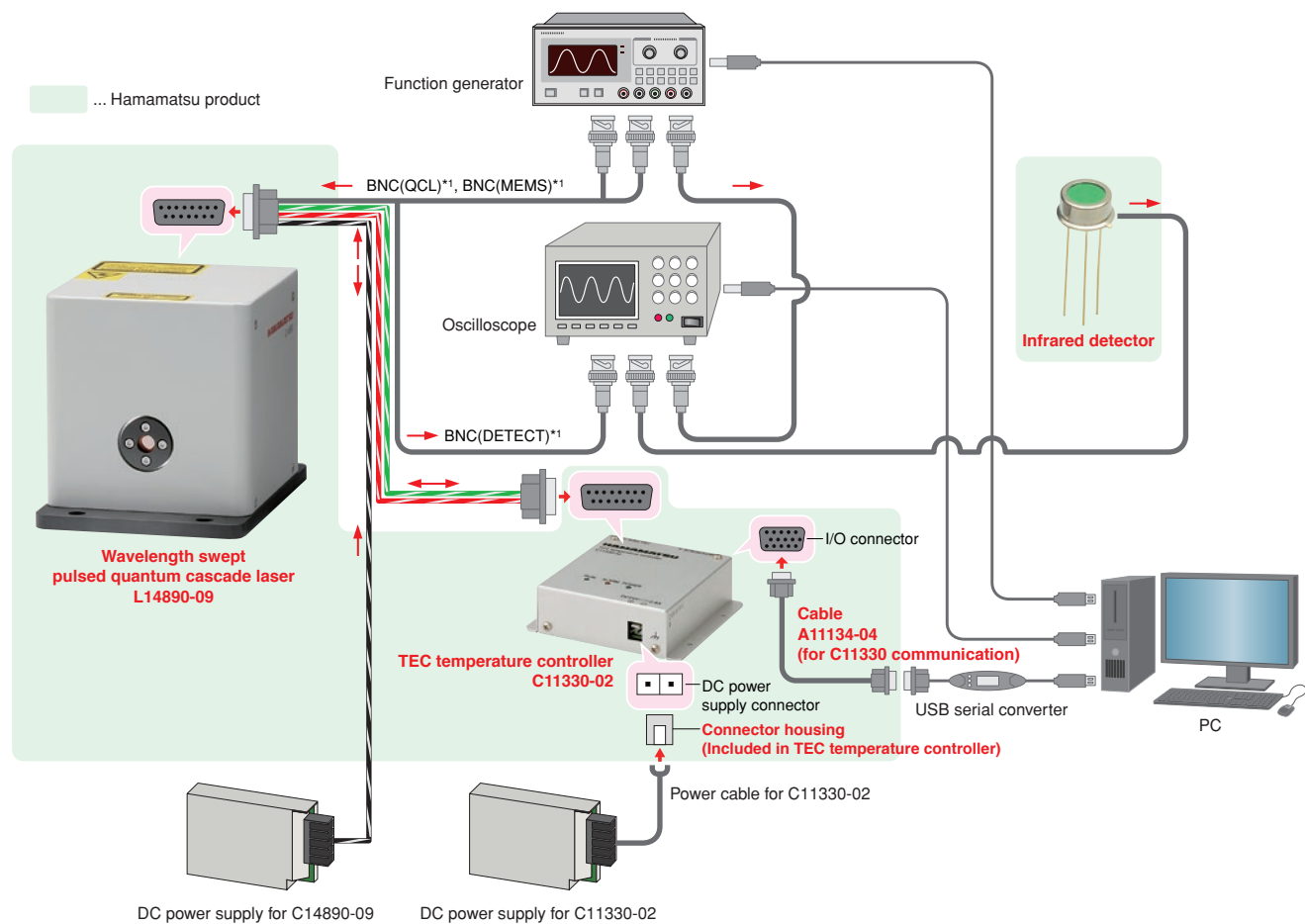

\*1: The name printed on the power cable attached to the product. Refer to "Power connector specifications" on page 4 for the functions.

\* The cable connected to the power connector is attached with the product. The cables connected to the DC power supply and the TEC temperature controller have to be manipulated by the user.

\* This configuration is recommended for working the sample software properly. Refer to the table of the required equipment for the model number of the recommended equipment.

### ● Accessories included product

Content	Quantity
Power cable (2 m)	1
Sample software (CD-R)	1
Instruction manual	1

### ● Necessary items for normal operation excluded product

Equipment	Manufacturer	Type No.
2-ch function generator	NF Corporation	WF1974
Oscilloscope	Keysight Technologies	DSOX2024A
DC Power supply	TDK lambda	HWS30A-24/A
TEC temperature controller	Hamamatsu Photonics K.K.	C11330-02 *1*2
PC	—	Windows 10 (64-bit) USB 2.0

\*1: A PC and an external DC power supply (24 V) are needed to control.


\*2: Recommended DC Power Supply: TDK lambda HWS50A-24/A

## ■ Thermistor

Characteristic	Symbol	Condition	Value	Unit
R25 resistance	R25	25 °C	10 ± 2.5 %	kΩ
Beta value	B	0 °C / 100 °C	3450	K

Wavelength swept pulsed quantum cascade laser L14890-09

Dimensional outline (unit: mm)

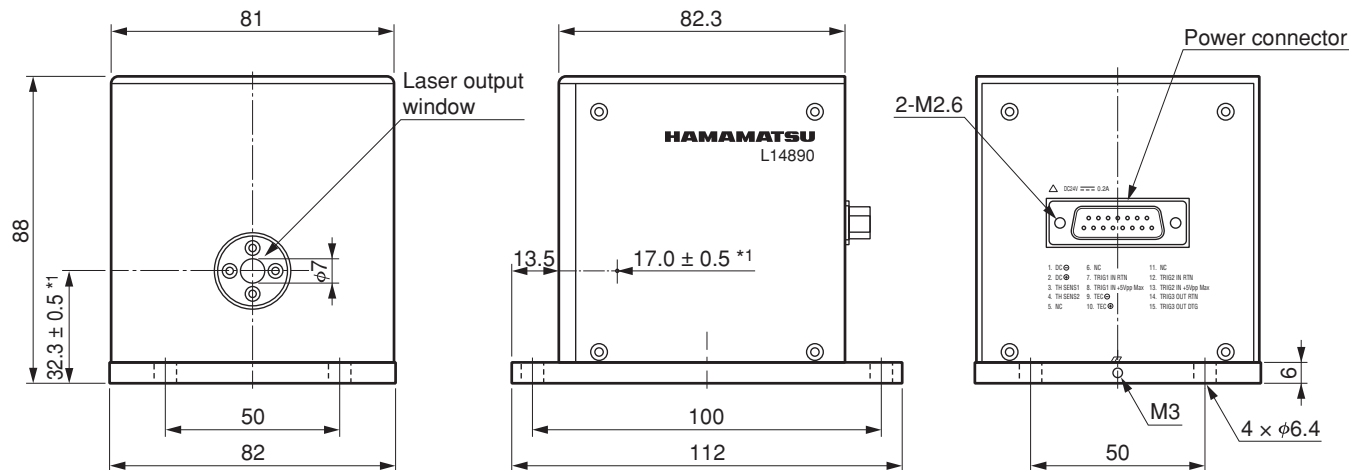


\*1: Emitter  
\* Tolerance: ±0.3 mm unless specified

LHJ3F0043-27\_A

Power connector specifications

15-pin D-sub connector (male)

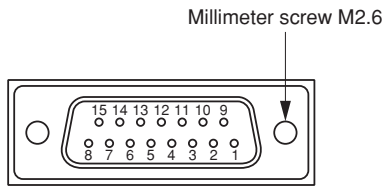


Power connector			Power cable
Pin No.	Symbol	Function	Wire color/name
1	24 V (-)	External power input	Black
2	24 V (+)	External power input	Black/White
3	Thermistor A	Temperature sensor output	Red
4	Thermistor B	Temperature sensor output	Red/White
5	N.C.	—	—
6	N.C.	—	—
7	QCL trig (-)	QCL operation signal input (Rtn)	BNC (QCL)
8	QCL trig (+)	QCL operation signal input (Sig)	
9	Peltier (-)	Peltier power input *1	Green
10	Peltier (+)	Peltier power input *1	Green/White
11	N.C.	—	—
12	MEMS (-)	MEMS operation signal input(Rtn)	BNC (MEMS)
13	MEMS (+)	MEMS operation signal input (Sig)	
14	DETECT (-)	MEMS detection signal input (Rtn)	BNC (DETECT)
15	DETECT (+)	MEMS detection signal input (Sig)	

\*1: The forward direction shows the cooling mode.  
\* The cable connected to the power connector is attached with the product.  
The cables connected to the DC power supply and the TEC temperature controller have to be manipulated by the user.

CLASS 3B LASER  
Invisible Laser Radiation: Avoid Exposure to Beam

•The Laser emits invisible laser radiation.  
The instrument which used the LASER, operated under ordinary conditions, is classified as Class 3B according to the laser product classification code IEC 60825-1. See IEC 60825-1, -14 for more details and safety operation concerning the above countermeasures.

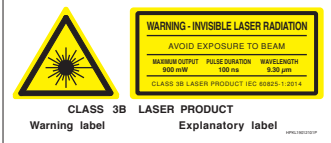

Information described in this material current as of October 2023. Specifications are subject to change without notice.

HAMAMATSU PHOTONICS K.K. www.hamamatsu.com

HAMAMATSU PHOTONICS K.K., Solid State Division  
1126-1 Ichino-cho, Higashi-ku, Hamamatsu City, 435-8558 Japan, Telephone: (81)53-434-3311, Fax: (81)53-434-5184

U.S.A.: HAMAMATSU CORPORATION: 360 Foothill Road, Bridgewater, NJ 08807, U.S.A., Telephone: (1)908-231-0960, Fax: (1)908-231-1218 E-mail: usa@hamamatsu.com  
Germany: HAMAMATSU PHOTONICS DEUTSCHLAND GMBH.: Arzbergerstr. 10, 82211 Herrsching am Ammersee, Germany, Telephone: (49)8152-375-0, Fax: (49)8152-265-8 E-mail: info@hamamatsu.de  
France: HAMAMATSU PHOTONICS FRANCE S.A.R.L.: 19 Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 10 E-mail: info@hamamatsu.fr  
United Kingdom: HAMAMATSU PHOTONICS UK LIMITED: 2 Howard Court, 10 Tewin Road, Welwyn Garden City, Hertfordshire, AL7 1BW, UK, Telephone: (44)1707-294888, Fax: (44)1707-325777 E-mail: info@hamamatsu.co.uk  
North Europe: HAMAMATSU PHOTONICS NORDEN AB: Torshamnsgatan 35, 16440 Kista, Sweden, Telephone: (46)8-509-031-00, Fax: (46)8-509-031-01 E-mail: info@hamamatsu.se  
Italy: HAMAMATSU PHOTONICS ITALIA S.R.L.: Strada della Moia, 1 int. 6 20044 Arese (Milano), Italy, Telephone: (39)02-93 58 17 33, Fax: (39)02-93 58 17 41 E-mail: info@hamamatsu.it  
China: HAMAMATSU PHOTONICS (CHINA) CO., LTD.: 1201, Tower B, Jiaming Center, 27 Dongsanhuan Beilu, Chaoyang District, 100020 Beijing, P.R. China, Telephone: (86)10-6586-6006, Fax: (86)10-6586-2866 E-mail: hpc@hamamatsu.com.cn  
Taiwan: HAMAMATSU PHOTONICS TAIWAN CO., LTD.: 13F-1, No.101, Section 2, Gongdao 5th Road, East Dist., Hsinchu City, 300046, Taiwan(R.O.C) Telephone: (886)3-659-0080, Fax: (886)3-659-0081 E-mail: info@hamamatsu.com.tw

Cat. No. LQCL3002E06  
OCT. 2023 IP