

## 38mm(1-1/2 Inch) Diameter, Extended Red Multialkali Photocathode Spectral Response 300 to 900 nm

### GENERAL

Parameter		Description/Value	Unit
Spectral Response		300 to 900	nm
Wavelength of Maximum Response		600	nm
Photocathode	Material	Extended Red Multialkali	–
	Minimum Effective Area	34	mm dia.
Window Material		Borosilicate	–
Dynode	Structure	Circular-cage	–
	Number of Stages	10	–
Direct Interelectrode Capacitances	Anode to Last Dynode	3	pF
	Anode to All Other Electrodes	4	pF
Suitable Socket		E678-12A(supplied)	–

### MAXIMUM RATINGS (Absolute Maximum Values)

Parameter		Value	Unit
Supply Voltage	Between Anode and Cathode	1500	Vdc
	Between Anode and Last Dynode	250	Vdc
Average Anode Current		0.2	mA
Ambient Temperature		-30 to 50	°C

### CHARACTERISTICS (at 25°C)

Parameter		Min.	Typ.	Max.	Unit
Cathode Sensitivity	Luminous (2856K)	120	200	–	μA/lm
	Red/White Ratio (R-68 filter)	–	0.3	–	–
Anode Sensitivity	Luminous (2856K)	20	50	–	A/lm
Gain		–	$2.5 \times 10^5$	–	–
Anode Dark Current (after 30min. storage in darkness)		–	8	30	nA
Time Response	Anode Pulse Rise Time	–	2.8	–	ns
	Electron Transit Time	–	40	–	ns

**NOTE:** Anode characteristics are measured with the voltage distribution ratio shown below.

### VOLTAGE DISTRIBUTION RATIO AND SUPPLY VOLTAGE

Electrodes	K	Dy1	Dy2	Dy3	Dy4	Dy5	Dy6	Dy7	Dy8	Dy9	Dy10	P
Ratio	2	1	1	1	1	1	1	1	1	1	1	1

Supply Voltage : 1000Vdc, K : Cathode, Dy : Dynode, P : Anode

# PHOTOMULTIPLIER TUBE R2066

Figure 1: Typical Spectral Response

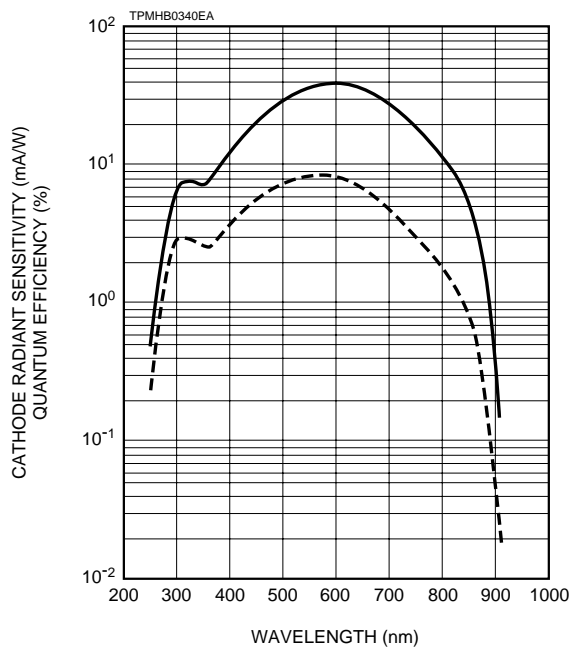


Figure 2: Typical Gain and Anode Dark Current


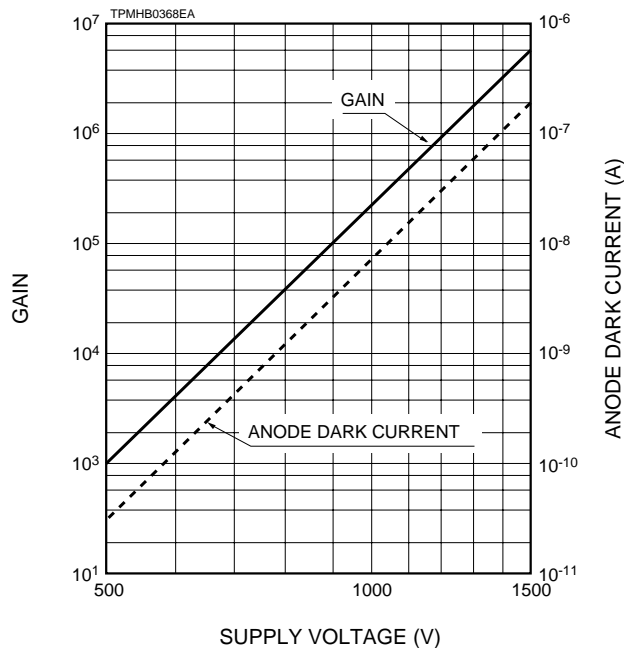


Figure 3: Dimensional Outline and Basing Diagram (Unit: mm)

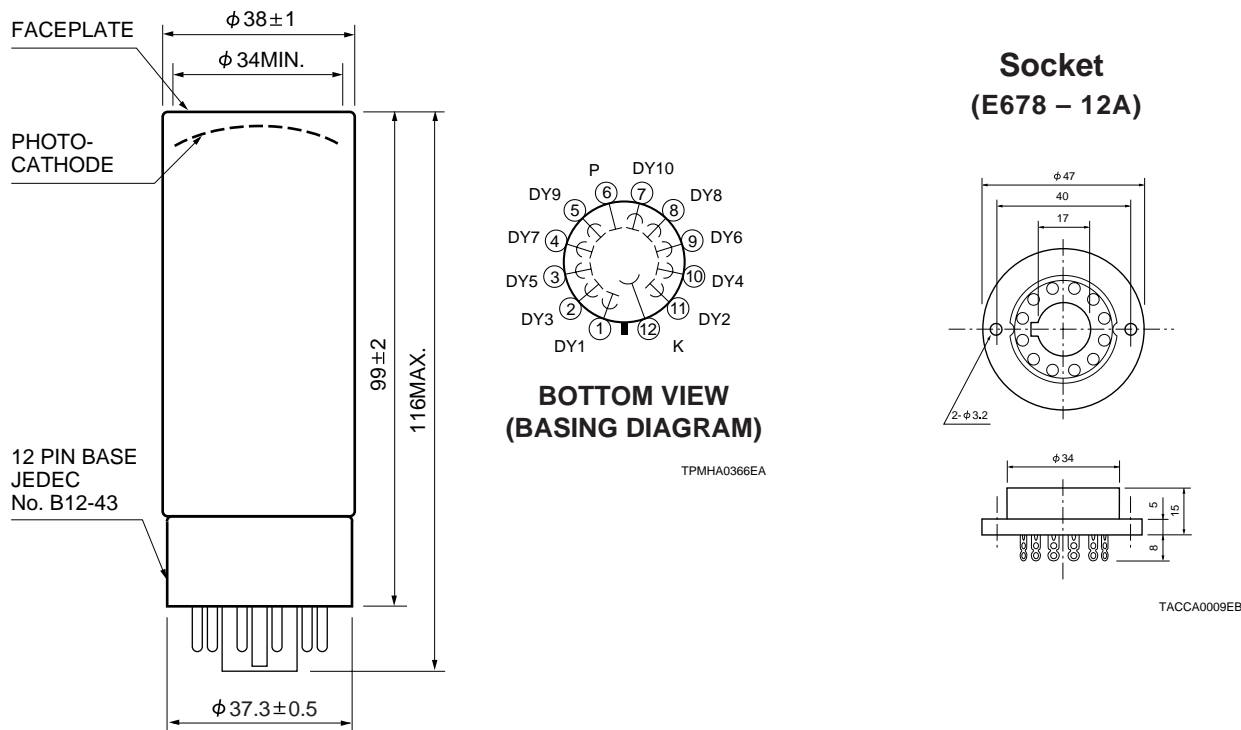

# HAMAMATSU

HOME PAGE URL <http://www.hamamatsu.com>

HAMAMATSU PHOTONICS K.K., Electron Tube Center

314-5, Shimokanzo, Toyooka-village, Iwata-gun, Shizuoka-ken, 438-0193, Japan, Telephone: (81)539/62-5248, Fax: (81)539/62-2205

U.S.A.: Hamamatsu Corporation: 360 Foothill Road, P. O. Box 6910, Bridgewater, N.J. 08807-0910, U.S.A., Telephone: (1)908-231-0960, Fax: (1)908-231-1218

Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49)8152-375-0, Fax: (49)8152-2658

France: Hamamatsu Photonics France S.A.R.L.: 8, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 10

United Kingdom: Hamamatsu Photonics UK Limited: Lough Point, 2 Gladbeck Way, Windmill Hill, Enfield, Middlesex EN2 7JA, United Kingdom, Telephone: 44(20)8-367-3560, Fax: 44(20)8-367-6384

North Europe: Hamamatsu Photonics Norden AB: Smidesvägen 12, SE-171-41 SOLNA, Sweden, Telephone: (46)8-509-031-00, Fax: (46)8-509-031-01

Italy: Hamamatsu Photonics Italia: S.R.L.: Strada della Moia, 1/E, 20020 Arese, (Milano), Italy, Telephone: (39)02-935 81 733, Fax: (39)02-935 81 741

TPMH1148E02  
DEC. 1999