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Poropare™

取扱説明書 / User Manual Ver. 3.1

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Table of contents (English edition)

1 Notes on using this product.....	9
2 Safety precautions.....	10
3 Handling precautions	10
4 Product description.....	11
5 Handling method.....	11
5.1 Unpacking and storage.....	11
5.2 Taking the transfer plate out of the case	12
5.3 Transferring the sample	12
5.4 Attaching to the dedicated adapter	13
6 Disposal.....	13
7 When you are in need of help	13

1 Precaution upon using this product

- Before using this product be sure to read the latest version of the user manual (this manual) and follow the safety precautions to ensure correct use.
 - This product is intended for use for time-of-flight mass spectrometers (TOFMS) equipped with an ion source utilizing the matrix-assisted laser desorption/ionization (MALDI) technique. Keep this product out of the reach of persons not familiar with this product.
 - To obtain detailed information on how to install this product in your mass spectrometer and how to use it after installation, please contact the dealer where you purchased this product or consult the mass spectrometer manufacturer.
 - Maintenance of peripheral equipment including mass spectrometers and replacement of consumable parts associated with the use of this product shall be performed by the customer on their own responsibility. Please note that mass spectrometry imaging measurements generally cause wear in mass spectrometers more rapidly than ordinary spectrometry.
 - This product was developed and manufactured to assist in sample preparation for mass spectrometry and its imaging. Do NOT use this product for any other applications.
 - This product is not guaranteed to be durable against any pretreatment.
 - Measurement results obtained using this product will vary depending on the method of use and the operating environment. This product is intended to assist in sample preparation and does not guarantee that the desired measurement results will be obtained.
 - We do not guarantee product's accuracy, completeness, merchantability, fitness for particular purpose or any other warranty, whether express or implied and we do not warrant that use of this product does not infringe intellectual property rights of third parties.
 - We cannot be held liable for any damages caused by (i) natural disasters such as earthquakes, lightning, windstorms and floods, (ii) fire (iii) enactment, amendment and repeal of laws, (iv) acts of third parties, (v) accidents during transport or storage, (vi) intentional or negligent misuse by the customer, (vii) usage under other unusual conditions, (viii) any other cases where we are not directly responsible for the damages.
 - To the maximum extent permitted under the applicable law, we assume no responsibility for any direct, indirect, consequential, special or incidental damages (including but not limited to failure of peripheral equipment including mass spectrometers, loss of business profits, business interruptions, loss of profits) related to or arising from the use or inability to use this product. In any given circumstances, our obligation to indemnify you for any damages resulting from or related to this product, either contractual, tort or otherwise, shall not exceed the amount equivalent to the price of this product.
 - Although every effort is made to ensure product quality, in the case of product damage caused during shipping or discoloration or deterioration (which affects measurement) found immediately after unpacking, we will replace the product with a new one, provided that all enclosed transfer plates are unused. In such a case, please contact the dealer where you purchased this product and then return the product along with the package described the lot number and the delivery slip or shipping information sheet. If no complaint is reported within 3 months from first receiving this product, the product will be deemed to have been accepted by the user as a good product.
-
- ✧ Reproduction or copying of this user manual in whole or in part is prohibited without our express permission.
 - ✧ The contents of this user manual are subject to change without prior notice.
 - ✧ The contents of this user manual are believed to be reliable. However, if you find any part of it unclear, then please contact us for assistance.

2 Safety precautions

Failure to observe the following precautions may result in injuries, equipment malfunctions or other unforeseen accidents.


- Do not use this product for samples that are generally considered hazardous.
Failure to do so may cause fire, burns, injuries, or equipment malfunctions. This product is not intended for use with samples that are explosive, ignitable, flammable, or otherwise hazardous.
- Do not apply excessive force to this product. Also, do not use this product under conditions where a localized force is applied (for example, due to using a pointed hard sample).
This product is made of glass so there is a risk of breakage.
- If this product is damaged, do not touch it with bare hands.
The broken surface is sharp and might cause injuries from cuts.
- Do not install this product in the equipment if damaged.
Debris might be left behind in the equipment causing it to malfunction.
- Always use an adapter (sold separately) dedicated for this product.
This product might otherwise drop off into the equipment, causing damage to the equipment.
- Securely install this product in place.
This product might otherwise drop off into the equipment, causing damage to the equipment.

3 Handling precautions


Failure to observe the following precautions may prevent this product from delivering full performance or might adversely affect the equipment.

- Avoid handling this product with bare hands and do not touch the active surface.
Unintended signals might be detected due to foreign matter or dirt that adheres to the product. When using gloves, use powder-free gloves.
- Do not reuse this product.
This product is disposable and cannot be reused. Reuse may generate unintended signals.
- Do not transfer highly adhesive samples or fragile, crumbly samples.
Excessive sample amounts adhering to the active surface of this product might prevent obtaining a signal. The vacuum in the mass spectrometer might also weaken.
- Store this product in a low-humidity environment and do not expose it to normal atmospheric conditions for long periods of time.
The active surface might lose its function or cause unintended signals to appear.
- Make sure there is electrical continuity between this product and the dedicated adapter.
Losing continuity between the product and dedicated adapter may cause failure to detect the signal or cause detection of unintended changes in the signal.

4 Product description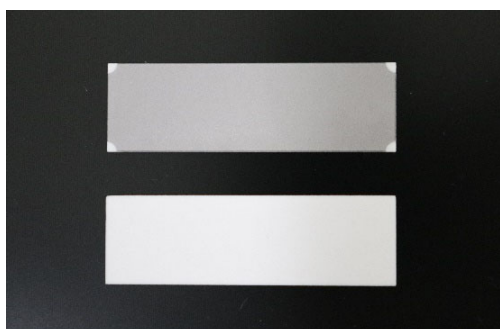

Top: Active surface (gray)
Bottom: Back side (white)

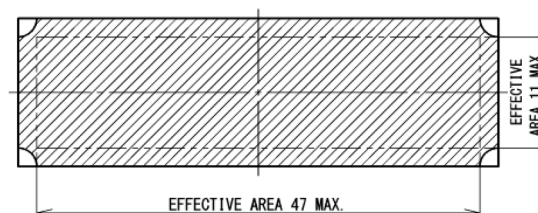Active area detailed view


Materials used	Soda glass, platinum (active surface)
Approximate product dimensions	52 mm on long side, 16 mm on short side, 1.05 mm (max.) in plate thickness
Maximum available active area	47 mm on long side, 11 mm on short side
Usage environment	Room temperature range (+10 °C to +30 °C)

5 Handling method

5.1 Unpacking and storage

Packaged state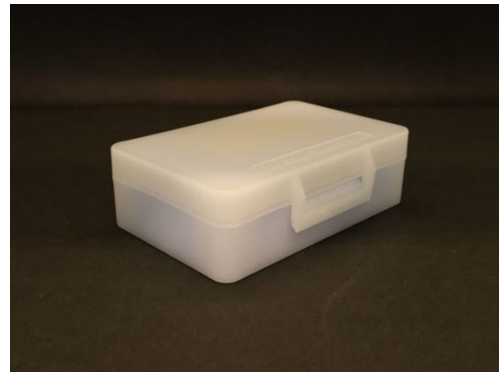


Case enclosed in package

When unpacking, be careful not to drop the case inside the package. This packaging is moisture-proofed to keep the product in a suitable environment during shipping. Please use the product as soon as possible after delivery. Be aware that once the package is opened, the moisture-proof effect will be lost, so open the package just before use. Since the case has no moisture-proofing, do not leave it in a room environment after taking it out of the package, and be sure to perform the sample transfer as soon as possible. When storing this product after unpacking, we recommend storing it in a hermetically sealed environment filled with a low-moisture inert gas such as in a dehumidified nitrogen desiccator. However, this is not a guarantee of product performance after unpacking and storing the product.

5.2 Taking the transfer plate out of the case

Avoid handling the transfer plate including the edges with bare hands and take care not to touch the active surface. When taking a transfer plate out of the case, first check which side is the active surface and then, while holding the transfer plate by the edges with fingers, take it out from the slits holding it. At this point, be careful to avoid applying excessive force to the transfer plate.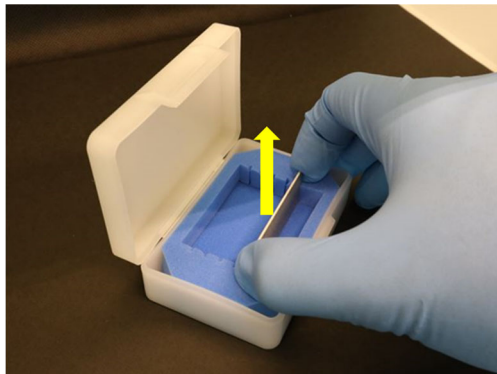


Wear powder-free gloves and take out the transfer plate by pulling straight up from the slits while being careful not to touch the active surface.


5.3 Transferring the sample

When transferring a sample, always comply with the instructions in "2. Safety precautions" and "3. Handling precautions." To prevent the transfer plate from bending or warping, place it on a flat hard surface so that the entire surface of the transfer plate is in uniform contact with the flat hard surface. The transfer method described in this section is just an example so it is not necessary to do exactly as described. However, please be careful to avoid cracking or damaging the transfer plate during the sample transfer. After transfer, check that there are no dust particles or foreign matter on the surface, and remove them if found.


Transfer process through contact with sample (Sample: Stains on film)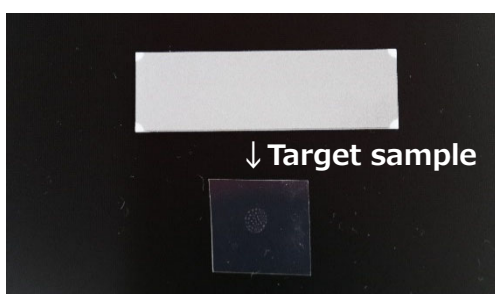


- (1) Place the transfer plate and sample on a flat hard surface.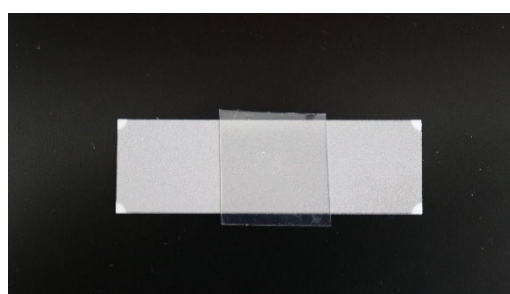

- (2) Make the active surface of the transfer plate comes in contact with the sample.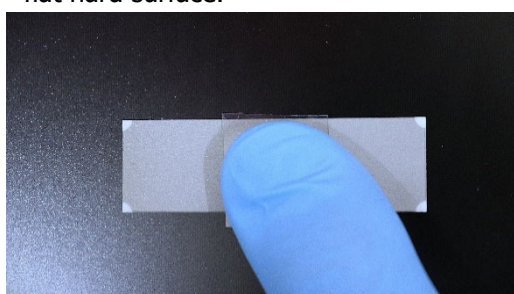

- (3) Press down on the transfer plate with your fingertip to make close contact with the sample.

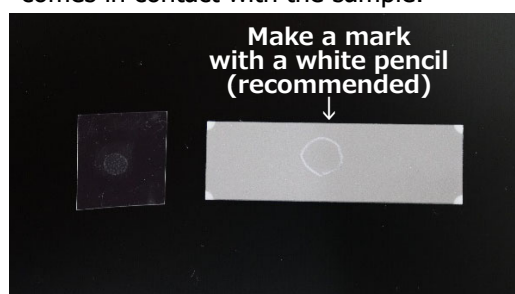- (4) Separate the transfer plate from the sample. (Marking the transfer point is recommended)

5.4 Attaching to the dedicated adapter

A dedicated adapter (sold separately) for Poropare™ transfer plates is required to install a transfer plate in a mass spectrometer. To find more information on the dedicated adapter, please contact the dealer where you purchased this product or consult the manufacturer of the mass spectrometer you are using. Securely attach the transfer plate to the dedicated adapter and make sure the active surface is aligned with the measurement surface. Also make sure there is electrical continuity between the active surface of the transfer plate and the dedicated adapter by using electrically conductive seals or other means. To find information on how to install the transfer plate in the mass spectrometer, follow the instructions in the mass spectrometer user manual. Do not fix the transfer plate on commercially available target plates or glass slides. The thickness of such plates or glass slides causes the measurement point to shift upward from its correct surface, increasing the risk of equipment failure.

After attaching the transfer plate to the dedicated adapter, be sure to remove any foreign matter and dust particles by using an air duster or similar device. Also be sure to check the transfer plate for any damage before installing it in the mass spectrometer. Set the measurement conditions according to the instructions in the mass spectrometer user manual. When adjusting the laser power, increase it gradually from the minimum output to a level that gives the desired signal. Also be careful not to make measurements with more laser power and number of laser shots than actually necessary.

6 Disposal

When disposing of this product, take appropriate measures in compliance with applicable regulations regarding waste disposal and correctly dispose of it yourself or entrust proper disposal to a licensed industrial waste disposal company. In any case, be sure to comply with the regulations in your country or state to ensure correct disposal.

7 When you are in need of help

To find more information on Poropare™, please access our website.

(<https://www.hamamatsu.com/all/en/product/optical-components/transfer-plate.html>)

There you can see a list of Q&A and refer to it as needed.

To find more detailed information on mass spectrometers, please contact the equipment manufacturer.

Reproduction or copying of this user manual in whole or in part is prohibited without our express permission.

The contents of this user manual are believed to be reliable. However, if you find any part of it unclear, then please contact us for assistance.

The contents of this user manual are subject to change without prior notice.

HAMAMATSU PHOTONICS K.K. www.hamamatsu.com

Electron Tube Division

314-5, Shimokanzo, Iwata City, Shizuoka Pref., 438-0193, Japan, Telephone: (81)539/62-5248, Fax: (81)539/62-2205

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