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INSTRUCTION MANUAL

C17198 series

- Before installing and using this product, be sure to read the “Safety Precautions”.
- This product undergoes complete in-house testing before shipment. However, unpack it carefully and check that all components and accessories are included and that there is no damage on the exterior or operational failure. If any components are missing or operational failure is suspected, contact your Hamamatsu sales office.

HAMAMATSU
PHOTON IS OUR BUSINESS

Table of Contents

Table of Contents	17
1. Before Using This Product	18
1.1 Safety precautions (Always ensure safety!)	18
1.2 Label positions	21
1.3 Features.....	21
1.4 Part names and accessories.....	22
2. Connections	24
Power connections	25
3. How to Use the C17198 series	26
3.1 Checking the connections	26
3.2 Using with Power supply	26
4. Precaution.....	27
4.1 Be sure to observe this precaution	27
4.2 Handling	27
4.3 Warranty	27
4.4 Aftersales service	27
5. Specifications.....	28

1. Before Using This Product

1.1 Safety precautions (Always ensure safety!)

This product was designed with safety and reliability in mind. However, like most electrical products, improper handling or mis operation may cause fire or electrical shock, resulting in accidents that may include serious injury or death. Be sure to observe the precautions listed in this manual or on the warning labels to prevent accidents.

Observe caution items for safety

Read the caution items described in this manual carefully and follow the instructions.

These caution items include general precautions relating to this product.

If any abnormality occurs ...

For example, if smoke, an unusual sound or odor is sensed, immediately turn off the power supply.

- Warning information in this manual is shown classified into the following items and by the extent of danger or damage that may result if the product is improperly used.

	WARNING	Failure to follow WARNING instructions could result in serious injury or death to the operator or person servicing the product
	CAUTION	Failure to follow CAUTION instructions may result in injury to the operator or the person servicing product, or damage to the product or peripheral equipment.

- Warning symbols used in this manual are classified as explained below. Make sure that you thoroughly understand the meaning of each symbol and follow the instruction. (These symbols are just examples.)

 CAUTION	Symbols showing a caution or warning you must pay attention to
 Must NOT do	Symbols showing what you must NOT do
 Must DO	Symbols showing what you must DO
	This symbol shows DC.

If the equipment is used in a manner not specified by this instruction manual, the protection provided by the equipment may be impaired. We are not liable for damage or losses resulting from failure to comply with precautions or instructions listed in this manual.

WARNING

Do not allow water or foreign matter to enter the product.

Fire

Electrical Shock

Must NOT do

Avoid placing any containers holding water near the equipment.

If water spills and enters the interior of the equipment, immediately turn off the power supply, then contact your Hamamatsu sales office correct servicing.

Do not disassemble or alter any part of this product.

Fire

Electrical Shock

Do NOT disassemble

Do not open the case of this product. A high voltage is applied to some internal parts of this product. Touching them may cause fire, electrical shocks or injury.

If smoke, an unusual odor or noise is coming from the product, immediately disconnect the power cable plug.

Fire

Electrical Shock

Must DO

Continuous operation may generate heat due to electrical shorts or insulation failure, resulting in fire or electrical shocks. Immediately turn off the power switch and disconnect the power cable plug.

Stop using the cables if damaged and disconnect the plugs.

Fire

Electrical Shock

Must DO

Continuous operation may generate heat due to electrical shorts or insulation failure, resulting in fire or electrical shocks. Immediately turn off the power switch and disconnect the power cable plug.

Remove any dust on the power cable plug.

Fire

Electrical Shock

Must DO

Dust or metallic chips adhering to the power cable plug may cause due to electrical shorts or insulation failure by moisture, resulting in fire or electrical shocks. If they are deposited, disconnect the plug and wipe off with dry cloth. Check to see if dust and foreign matter adhere to the plug occasionally.

Follow the authorized disposal methods if discarding this product.

Must DO

Please follow the applicable regulations regarding disposal of hazardous materials and industrial wastes in your country, state, region or province.

Do not touch the power cable plug with wet hands.

Electrical Shock

Must NOT do

Electricity conducts through water so touching the power cable plug with wet hands may result in electrical shocks.

CAUTION

Avoid installing this product in locations subject to excessive soot, steam, humidity, dust and vibration.

Fire

Must NOT do

Soot, steam, humidity, dust, corrosive gases and inflammable gases may harm this product. Especially, if dust adheres to the input window of this product, transmittance may degrade significantly.

Do not tug on the power cable to unplug it.

Fire

Electrical Shock

Must NOT do

Tugging on the power cable may damage the cable. Continuous operation with a damaged cable may generate heat by due to electrical shorts or insulation failure, resulting in fire or electrical shocks.

Unplug the power cable for safety during maintenance or long-term storage.

Fire

Electrical Shock

Must DO

Touching an internal part by mistake may result in electrical shocks. Storing the product left energized for extended periods of time may cause fire.

Read this manual thoroughly and follow the instructions.

Fire

Electrical Shock

Must DO

If the equipment is used in a manner not specified by this instruction manual, the protection provided by the equipment may be impaired. We are not liable for damage or losses resulting from failure to comply with precautions or instructions listed in this manual.

Use a relay ring that matches the C-mount component to be attached to the input window.

Must DO

If you want to attach a C-mount component which requires a threaded portion longer than 5 mm, then use a commercial C-mount relay ring having a correct length. If you attempt to forcefully attach a C-mount component which does not match the threaded portion, that component or the electromagnetic wave input window of the C17198 series might be damaged. Always use a proper C-mount adapter.

Avoid using this product in a vacuum.

Fire

Electrical Shock

Must NOT do

If this product is used in a vacuum, an internal component at a high voltage may cause discharge, resulting in fire or electrical shocks.

Disconnecting device.

Fire

The device corresponding to the disconnecting device (power shutoff device) for this product is the AC Adapter (that connects to the power outlet). In case of an emergency that requires the power to be shut off immediately to avoid danger, unplug the AC Adapter from the power outlet. When installing this product, ensure enough space so that the AC Adapter can easily be unplugged in case of emergency.

1.2 Label positions

Power supply / C17371

C17198 series

Install this product so that warning labels affixed on the unit can be easily seen during operation. If the label peels off, becomes dirty, or illegible, replace it with a new label. For a new warning label (available at an extra charge), Please contact our sales office or sales representative.

1.3 Features


The C17198 series is a THz-wave sensor module incorporating a THz Image Intensifier.

<Applications>


- Real time THz-wave beam monitor
- THz transmission imaging

1.4 Part names and accessories

When the product is delivered to you, unpack it carefully and check that there is no damage on the exterior and that all components and accessories are included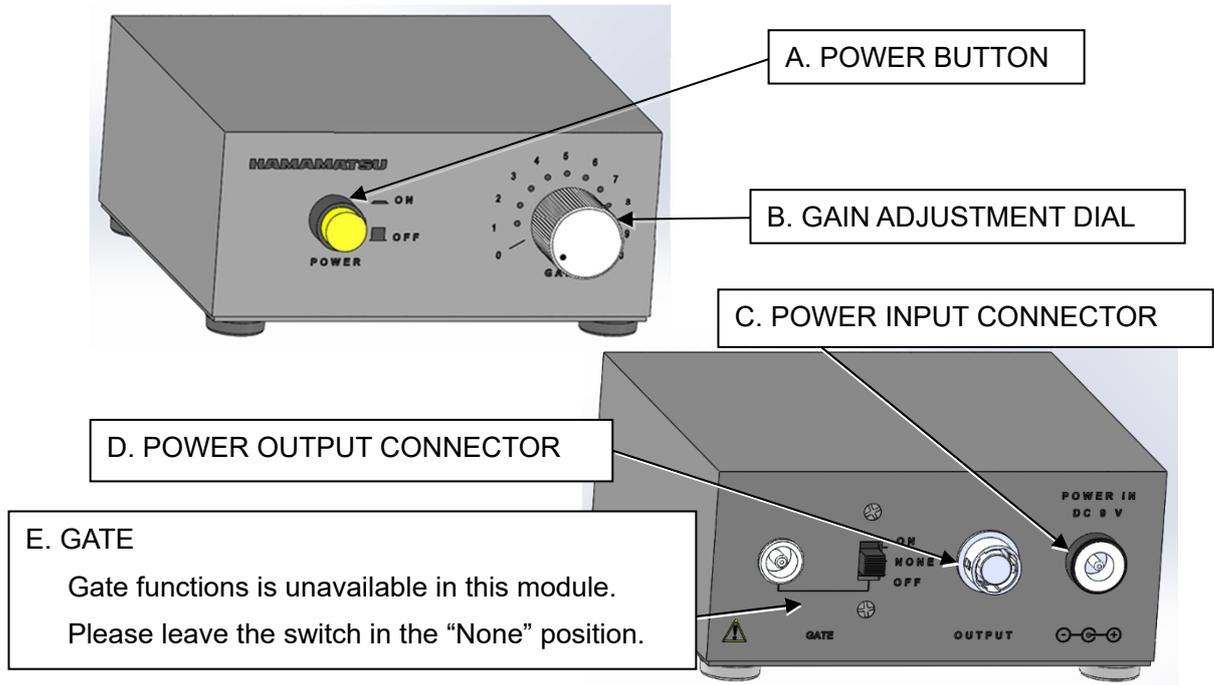

Power supply/C17371

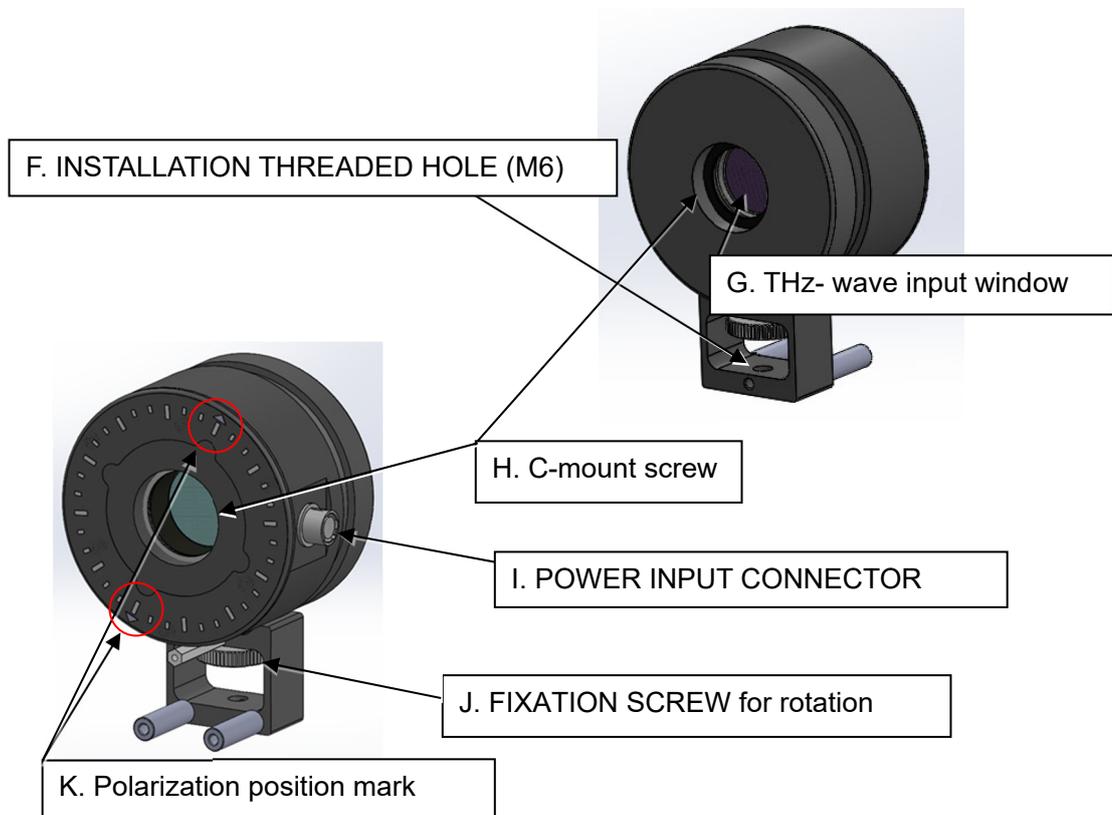C17198 series

A. POWER BUTTON

When the power is on, the LED lights up.

When a battery pack is used, the LED light intensity will decrease and the LED will blink when the voltage drops.

If the light intensity drops or blinks, stop using the battery pack and replace it with a new battery.

B. GAIN ADJUSTMENT DIAL

In case that the gain control dial is 0, the gain is minimum.

C. POWER INPUT CONNECTOR

This is for power input to the power supply.

D. POWER OUTPUT CONNECTOR (6-pin connector)

This is for the power output of the C17198 series.

E. GATE

Gate functions is unavailable in this module. Please leave the switch in the "None" position.

F. INSTALLATION THREADED HOLE (M6)

The standard of the screw is M6. Use this hole for mounting optical system on the optical table.

G. THz wave input window

THz-wave enters the sensitive area through this window.

H. C-mount screw

This is a standard screw-threaded mount and allows attaching various C-mount components such as a lens, diaphragm and shutter.

I. POWER INPUT CONNECTOR (6-pin connector)

This is the C17198 series for power input.

J. FIXATION SCREW for rotation

This is to unscrew it and can turn a device.

K. Polarization position mark

This mark indicates the polarization direction confirmed at the factory.


■Package contents

					
C17198 series 1 piece	Power supply/ C17371 1 piece	**1AC Adapter 1 piece	Power cable (120 cm) 1 piece	Battery box (9V battery) 1 piece	Safety Precautions 1 vol.

**1Not included in C17371-00.


2. Connections

However, some setups are required before beginning measurement. This section explains how to make setups for measurement in the following points.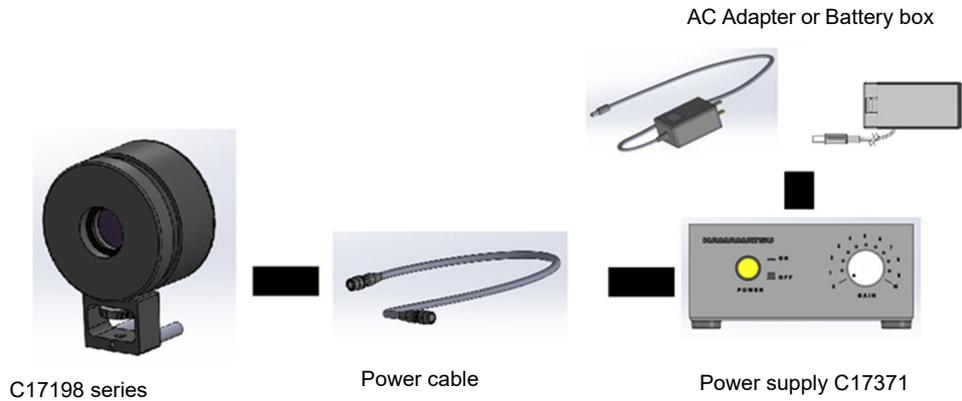

Connection block diagram at AC adapter

1. Connect the Power cable to the C17198 series and Power supply C17371.
2. Connect the AC adapter to the Power supply C17371.
3. Plug the AC Adapter into the power outlet.

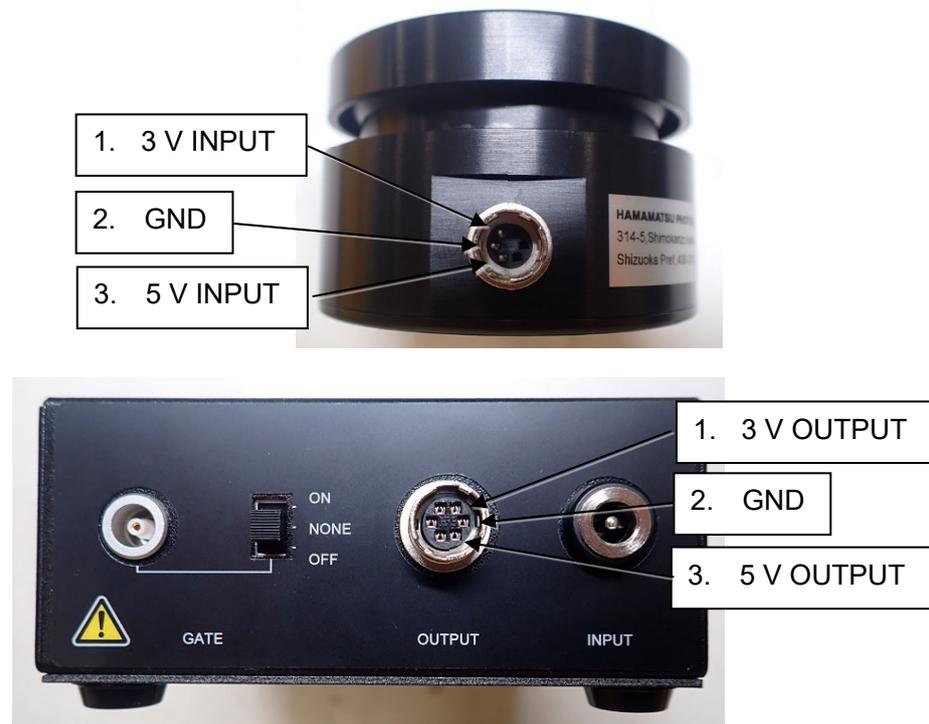
Connection block diagram at Battery box

1. Connect the Power cable to the C17198 series and Power supply C17371.
2. Open the battery box lid and connect a 9V dry cell battery to the socket inside.
3. Connect the Battery box to the Power supply C17371.
4. If the LED blinks or lights up but the light intensity is significantly reduced when the power button is pressed, discontinue use and replace the battery with a new one, as the battery is suspected to be depleted.

Battery box open

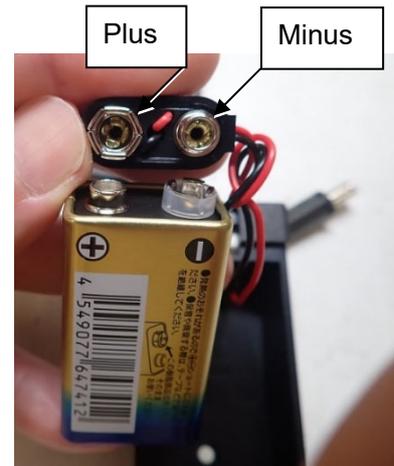Aspect of connection


CAUTION: This product is designed for indoor use and cannot be installed and used outdoors or in locations exposed to wind and rain.

WARNING: Use the AC Adapter and Power cable supplied with this product.

Power connections

The POWER connectors on the side of the C17198 series are used to supply power to the C17198 series. To connect to this POWER connectors, align the guide mark on the plug (supplied) with the POWER connectors, and insert it into position while holding the plug end. To disconnect, hold the ring on the plug and pull outwards.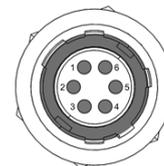

Example. HR10A-7R-6SC Plug

3. How to Use the C17198 series

3.1 Checking the connections

Before turning the power on, check the following points.

Connection of measurement systems:

Check the connections of the measurement systems. If the connections are incomplete and light leaks, accurate measurement can not be performed and these situations may even damage equipment.

3.2 Using with Power supply

Turning the power on or off

To turn the power on:


1. Check that the upper window of the control voltage adjustment dial reads "0". If not "0", turn the dial counterclockwise to return it to "0".
2. Turn on the POWER BUTTON of the Power supply.
The POWER BUTTON LED lights up and power is supplied to the C17198 series.
3. Turn the GAIN ADJUSTMENT DIAL clockwise to adjust the sensitivity of the C17198 series.

To turn the power off:

1. Turn the GAIN ADJUSTMENT DIAL counterclockwise to return it to "0".
2. Turn off the POWER BUTTON of power supply.
The POWER BUTTON LED also turns off.

3.3 Checking the Polarization direction

The sensitivity is maximized when the beam amplitude direction is aligned with the Polarization position mark.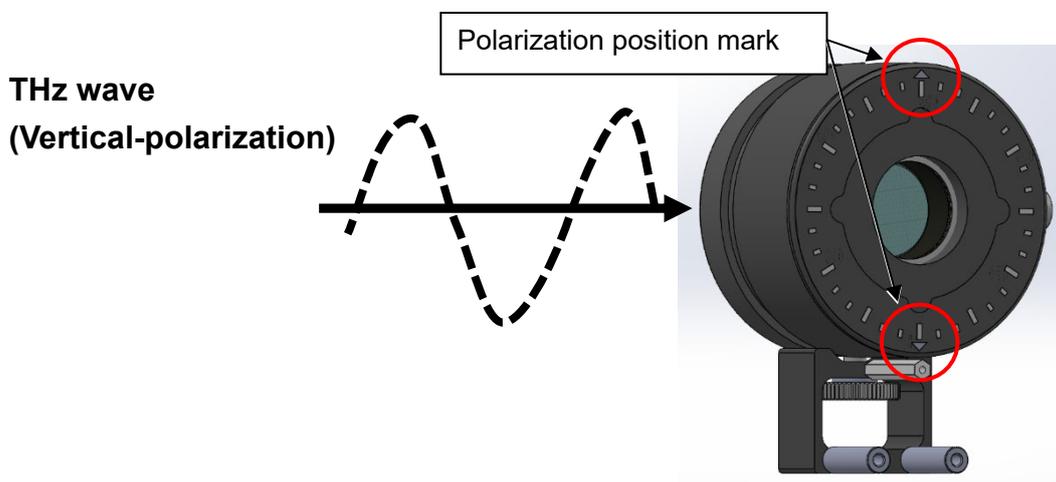

4. Precaution

4.1 Be sure to observe this precaution

When this product is not used for long periods of time, cover it with a plastic cover and store in a place with low humidity and no exposure to light.

4.2 Handling

Use sufficient caution to keep the input window clean. If dust or dirt gets on the window, use a soft cloth moistened with ethyl alcohol and gently wipe it off. Avoid using any organic solvents except ethyl alcohol.

4.3 Warranty

This product is warranted for one year from the date of delivery, whichever comes first. If any failure is found in the workmanship or materials within this warranty period, Hamamatsu will repair or replace the defective parts without charge. The warranty shall not apply to failure in the following cases.

- (2) Failure or trouble was caused by mis operation, mishandling, or not following the instructions and precautions described in this manual.
 - (3) The product has been modified electrically or mechanically by the customer.
 - (4) Failure was caused by accidents such as natural or man-made disasters.
- The warranty is limited to repair or replacement.
-

4.4 Aftersales service

If a failure has occurred after extended periods of operation, which was caused by breakage or wear of replaceable parts, please ship the product back to us for repair, replacement or adjustment. (This cost will be charged to you.)

5. Specifications

GENERAL

Parameters		Ratings	Units
Cathode	TYPE	Metamaterial	-
	ID	01	-
	Minimum Effective Diameter	16.0	mm
Input Window	Material	Silica Glass	-
	Thickness	5.5	mm
MCP (Micro Channel Plate)		2 Stages	-
Output Window	Material	Fiber Optics Plate	-
Phosphor Screen	Material	P43	-
	Minimum Effective Diameter	16.0	mm

RATINGS

Parameters		Min.	Typ.	Max.	Units	
POWER INPUT	*1AC adapter	Input Voltage	100	-	240	Vac
		Input Current	-	-	0.4	A
		Input Rated frequency	50 / 60			Hz
	BATTERY BOX	Input Voltage	7.3	9	9.5	Vdc
		Input Current	-	-	60	mA
Ambient Temperature	Storage	-20	-	+50	°C	
	Operating	+5	-	+50	°C	
Ambient Humidity	Storage	-	-	85	%RH	
	Operating	-	-	80	%RH	
Recommended spectral response		0.7 to 1.3			THz	
Threshold E-field			5	10	kV/cm	
Polarization dependence		Yes			-	
Dimensions	C17198 series	66(W) × 92(H) × 43(D)			mm	
Weight	C17198 series	Approx.310			g	

*1Not included in C17371-00.

OTHERS

Location for use		Indor
Altitude for use		2000 meters or less
Overvoltage category		II
Pollutin degree		2
EMC compliance	IEC	IEC 61326-1 Emission: Group 1, Class B, Immunity: table 2 ^{*1}
	FCC ^{*2}	FCC47 CFR Part 15 Subpart B, Class B
Safety standard compliance		IEC/EN 61010-1
IP code		IP 20
RoHS category		9


^{*1}Performance Level

Criterion A	Phosphor Brightness	Fluctuations in fluorescence intensity should be within $\pm 60\%$.
Criterion B	Phosphor Brightness	Return to a normal state automatically after the test.
Criterion C	Phosphor Brightness	Return to a normal state by operator's manual operation after the test.

^{*2}FCC Note

<p>This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.</p> <p>This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:</p> <ul style="list-style-type: none"> • Reorient or relocate the receiving antenna. • Increase the separation between the equipment and receiver. • Connect the equipment into an outlet on a circuit different from that to which the receiver is connected. <p>Consult the dealer or an experienced radio/TV technician for help.</p>
--


Dimensions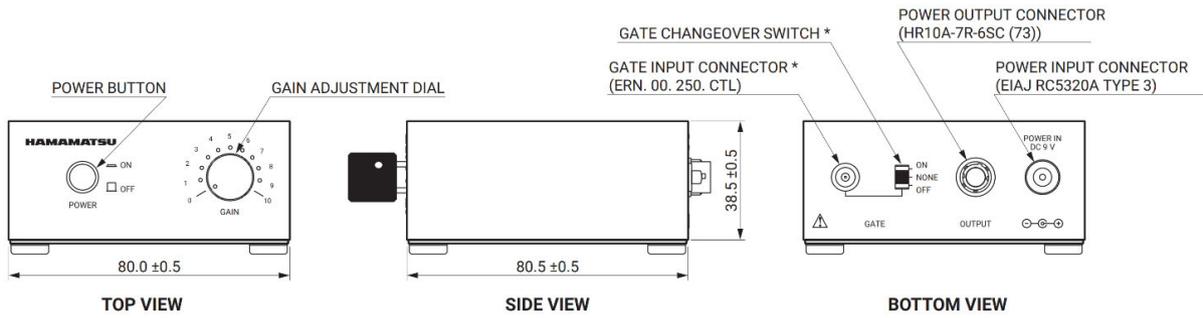

* Gate functions is unavailable in this module.

Power supply / C17371

Unit: mm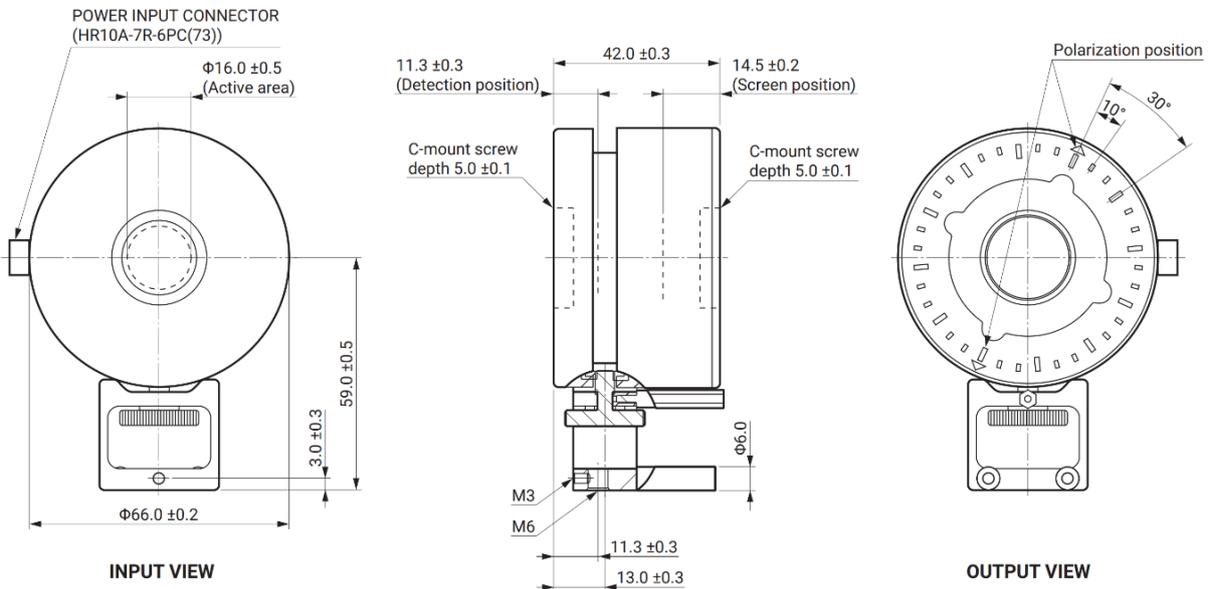

C17198 series

Unit: mm

@a1-0033-3 2025/05

HAMAMATSU PHOTONICS K.K. www.hamamatsu.com

Electron Tube Division

314-5, Shimokanzo, Iwata City, Shizuoka Pref., 438-0193, Japan, Telephone: (81)539/62-5248, Fax: (81)539/62-2205

U.S.A.: HAMAMATSU CORPORATION: 360 Foothill Road, Bridgewater, NJ 08807, U.S.A., Telephone: (1)908-231-0960, Fax: (1)908-231-1218

Germany: HAMAMATSU PHOTONICS DEUTSCHLAND GMBH: Arzbergerstr. 10, 82211 Herraching am Ammersee, Germany, Telephone: (49)8152-375-0, Fax: (49)8152-265-8 E-mail: info@hamamatsu.de

France: HAMAMATSU PHOTONICS FRANCE S.A.R.L.: 19 Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 00 E-mail: info@hamamatsu.fr

United Kingdom: HAMAMATSU PHOTONICS UK LIMITED: 2 Howard Court, 10 Tewin Road, Welwyn Garden City, Hertfordshire, AL7 1BW, UK, Telephone: (44)1707-294888, Fax: (44)1707-325777 E-mail: info@hamamatsu.co.uk

North Europe: HAMAMATSU PHOTONICS NORDEN AB: Torshamnsgatan 35, 16440 Kista, Sweden, Telephone: (46)8-509-031-00, Fax: (46)8-509-031-01 E-mail: info@hamamatsu.se

Italy: HAMAMATSU PHOTONICS ITALIA S.R.L.: Strada della Moia, 1 int. 6 20044 Arese (Milano), Italy, Telephone: (39)02-93 58 17 33, Fax: (39)02-93 58 17 41 E-mail: info@hamamatsu.it

China: HAMAMATSU PHOTONICS (CHINA) CO., LTD.: 1201, Tower B, Jiaming Center, 27 Dongsanhuang Bellu, Chaoyang District, 100020 Beijing, P.R. China, Telephone: (86)10-6586-6006, Fax: (86)10-6586-2866 E-mail: hpc@hamamatsu.com.cn

Taiwan: HAMAMATSU PHOTONICS TAIWAN CO., LTD.: 13F-1, No.101, Section 2, Gongdao 5th Road, East Dist., Hsinchu City, 300046, Taiwan (R.O.C) Telephone: (886)3-659-0080, Fax: (886)3-659-0081 E-mail: info@hamamatsu.com.tw