

Active Silicon Phoenix and FireBird (CameraLink)

DCAM Version

DCAM Module	19.12.322.5902	(for 32-bit)
	19.12.642.5902	(for 64-bit)
DRIVER	6.86.300.5902	(for PHX; D24CL-PE1)
	7.05.140.5902	(for FBD; 1xCLD-2PE8)
	8.13.300.5902	(for FBD; 2PE4, 2xCLD-2PE8)

Cards

Cards	Camera Link Configuration	PC Bus Type	Support OS	Note
AS-PHX-D24CL-PE1	Base	PCI Express x1	Windows 10 (*25) 32-bit / 64-bit (x64) Windows 7 / 8.1 32-bit / 64-bit (x64)	Half Length
AS-FBD-1XCLD-2PE8	Deca	PCI Express x8 Gen2 (*15)(*22)		Half Length / Dual Slot
AS-FBD-2XCLD-2PE8	Dual Deca			Half Length Low Profile PCB - Full Height Bracket
AS-FBD-1XCLD-2PE4L-F	Deca / Dual Base / Full / Medium	PCI Express x4 Gen2 (*15)		Half Length Low Profile PCB - Low Height Bracket
AS-FBD-1XCLD-2PE4L-L	Deca / Dual Base / Full / Medium	PCI Express x4 Gen2 (*15)		

Cameras

Fast speed CameraLink cameras

Cameras	Nickname	Supported Card	Note
C13440-20C(U)	ORCA-Flash4.0 (V3)	AS-FBD-1XCLD-2PE4L-F AS-FBD-1XCLD-2PE4L-L AS-FBD-2XCLD-2PE8	
C11440-22C(U)	ORCA-Flash4.0 (V2)	AS-FBD-1XCLD-2PE8 AS-FBD-2XCLD-2PE8	
C11440-10C	ORCA-Flash2.8	AS-PHX-D24CL-PE1	(*18)
C10000-A01	TDI Board Camera	AS-FBD-1XCLD-2PE4L-F AS-PHX-D24CL-PE1	
C10000-801	TDI Camera		

CameraLink cameras

Cameras	Nickname	Supported Card	Note
C8000-30		AS-PHX-D24CL-PE1	
C8484-xxC(P)			
C9100-13	ImagEM		
C9100-14	ImagEM 1K		
C12741-11	InGaAs VGA Camera	AS-FBD-1XCLD-2PE4L-F	(*New)
C9750-xxxx(N-C)	X-Ray Line	AS-PHX-D24CL-PE1	
C10400-xx			
C12450-27FGC-C			(*New)
C10650-xx	X-Ray TDI	AS-FBD-1XCLD-2PE4L-F	
C12200-321/461			
C12300-321			
C10800-xx-C	X-Ray Line Dual Energy	AS-PHX-D24CL-PE1	

Recommendation

It is highly recommended to disable C-state processor control in your PC's BIOS, else you may get sporadic corrupted images transferred to the PC. See Note (*23).

Active Silicon FireBird (CoaXPress)

DCAM Version

DCAM Module	19.12.642.5902	(for 64-bit)
DRIVER	8.26.700.5902	(for 4XCXP6-2PE8)
	8.26.500.5902	(for 2XCXP6-2PE8)

Cards

Cards	PC Bus Type	Support OS	Note
AS-FBD-4XCXP6-2PE8	PCI Express x8 Gen2	Windows 10 (*25) 64-bit (x64)	Half Length
AS-FBD-2XCXP6-2PE8	PCI Express x8 Gen2		Windows 7 / 8.1 64-bit (x64)

Cameras

CoaXPress cameras

Cameras	Nickname	Supported Card	Note
C14120-20P	ORCA-Lightning	AS-FBD-4XCXP6-2PE8	
C14440-20UP	ORCA-Fusion	AS-FBD-2XCXP6-2PE8	

Recommendation

It is highly recommended to disable C-state processor control in your PC's BIOS, else you may get sporadic corrupted images transferred to the PC. See Note (*23).

Active Silicon Phoenix (LVDS/RS422)

DCAM Version

DCAM Module	19.12.322.5902	(for 32-bit)
	19.12.642.5902	(for 64-bit)
DRIVER	6.86.400.5902	

Cards

Cards	PC Bus Type	Support OS	Note
AS-PHX-D36-PE1	PCI Express x1	Windows 10 (*25) 32-bit / 64-bit (x64) Windows 7 / 8.1 32-bit / 64-bit (x64)	Half Length

Cameras

LVDS/RS-422 cameras

Cameras	Nickname	Supported Cards	Note
C9750-xx	X-Ray Line	AS-PHX-D36-PE1	
C10400-xx			

Recommendation

It is highly recommended to disable C-state processor control in your PC's BIOS, else you may get sporadic corrupted images transferred to the PC. See Note (*23).

GigE

DCAM Version

DCAM Module	19.12.339.5902	(for 32-bit)
	19.12.659.5902	(for 64-bit)
DRIVER	5.1.104642.5902	

Cameras

Camera or Sensors	Nickname	Support OS	Note
C15333-10E	InGaAs Line Camera	Windows 10 (*25) 32-bit / 64-bit (x64) Windows 7 / 8.1 32-bit / 64-bit (x64)	(*New)
C12450-27FGC-G	X-Ray Line		(*New)
C12902D-40	Flat Panel Sensor		
C12903D-40			
C12504D-56			
C12505D-56			
C10500D-42/43/70			
C10502D-42/43/70			
C10900D-40			
C10901D-40			
C11700DK-40			
C11701DK-40			



Recommendations

See the next pages.

GigE – Cont'd

Recommendations


It is highly recommended to tweak these settings for the Gigabit Network Adapter for which the GigE device is connected, provided they are available for adjustment in Device Manager \ Properties for the adapter driver:

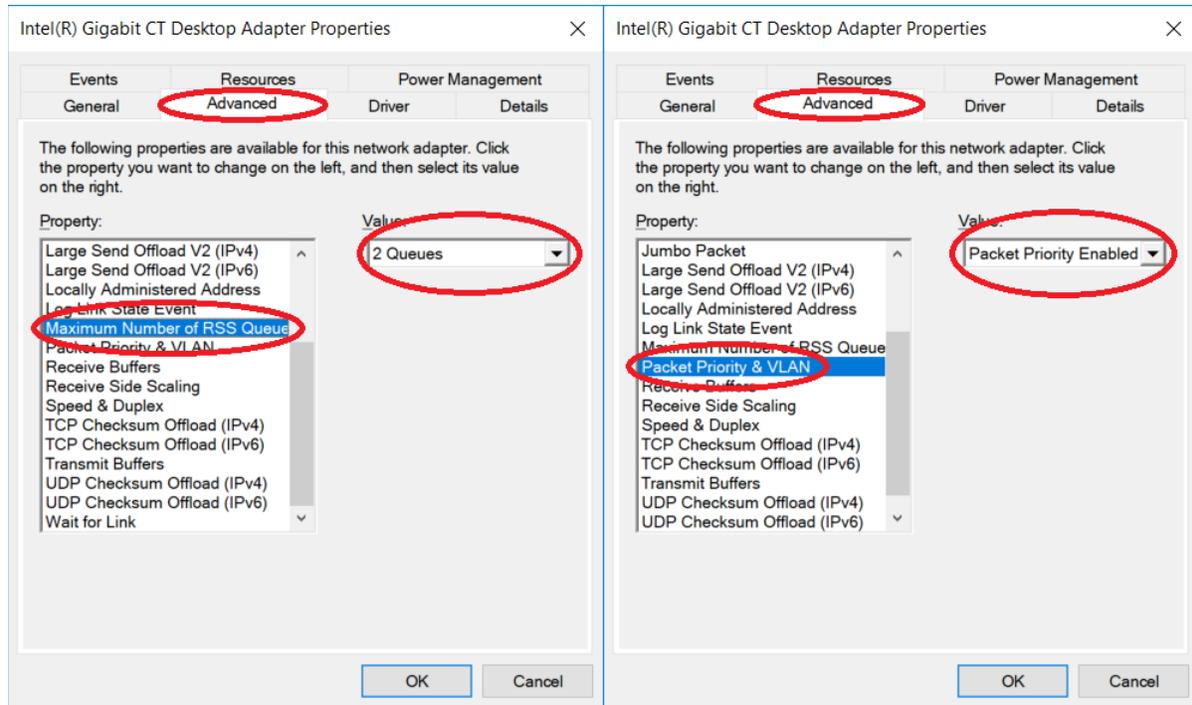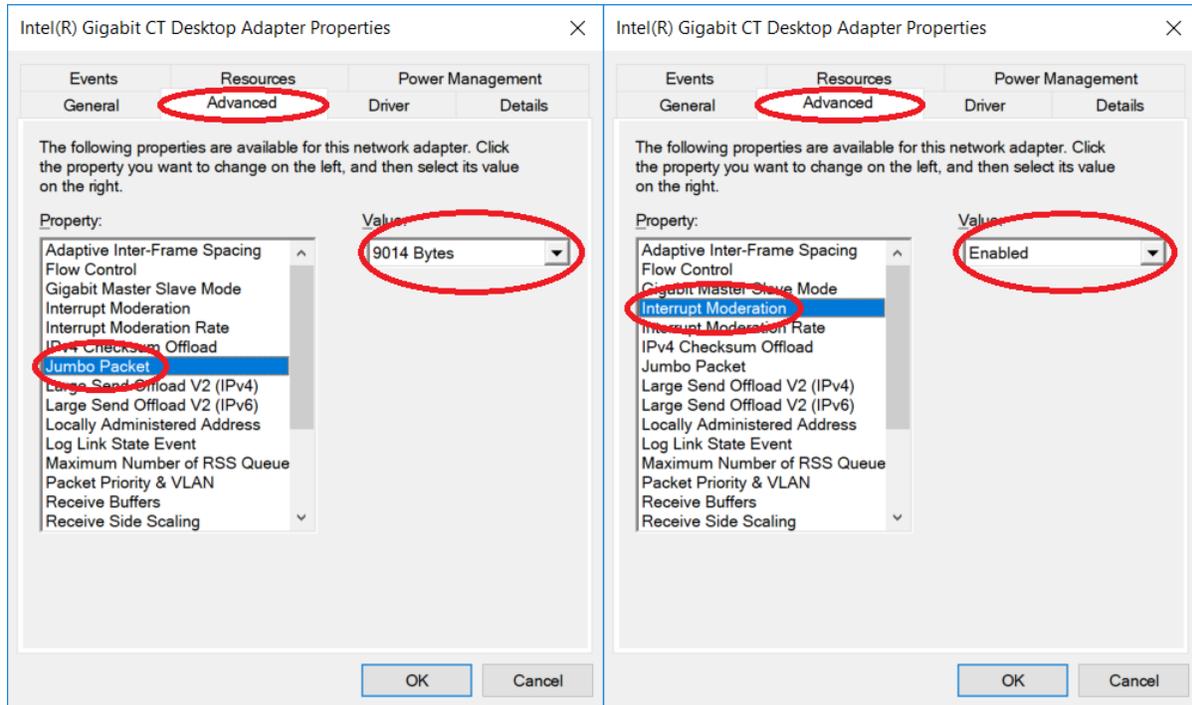

GigE – Cont'd

Recommendations

It is highly recommended to tweak these settings for the Gigabit Network Adapter for which the GigE device is connected, provided they are available for adjustment in Device Manager \ Properties for the adapter driver:

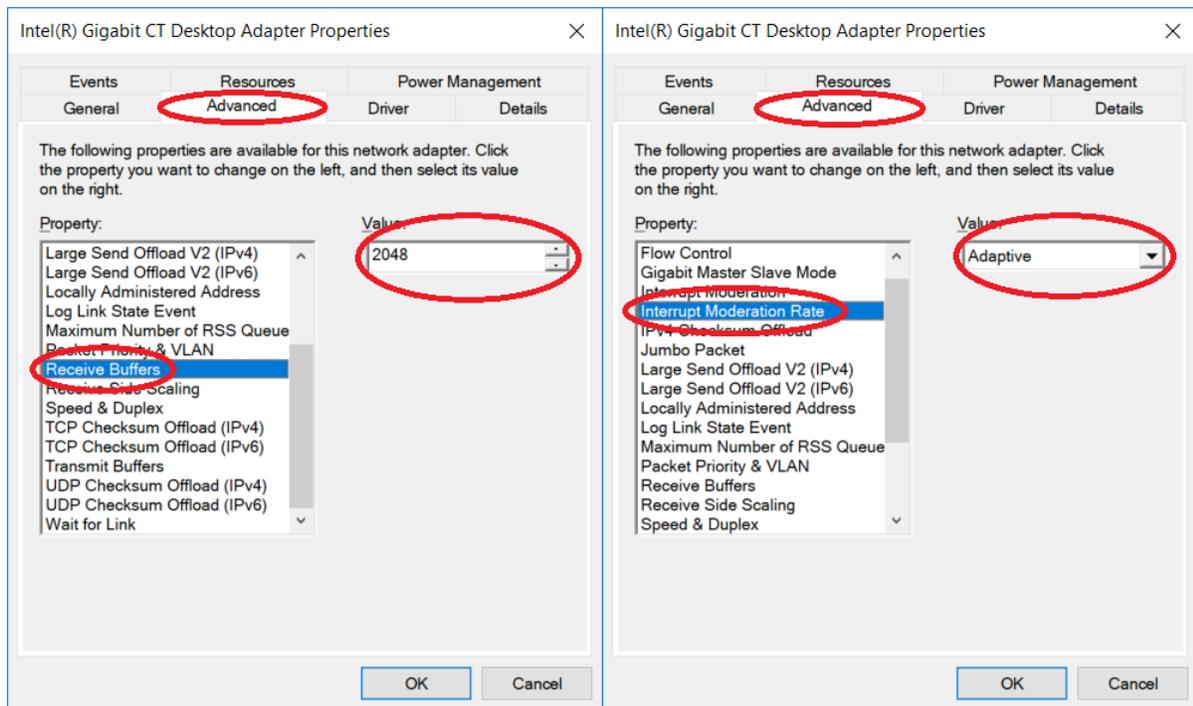

Notables:

- Jumbo Packets has the biggest effect to sustainable FPS and bandwidth. Set this setting to the highest possible by the adapter driver.
- Receive buffers should be set to maximum allowed by the driver if the setting exists.

USB

DCAM Version

DCAM Module	19.12.322.5902	(for 32-bit)
	19.12.642.5902	(for 64-bit)
DRIVER	1.2.6.5902	(for USB 3.0 and C10633)
	2.12.2.5902	(for others)

Cameras

Cameras	Nickname	USB2.0	USB3.0	Support OS	Note
C14440-20UP	ORCA-Fusion		✓	Windows 10 (*25) 32-bit / 64-bit (x64)	
C13440-20CU	ORCA-Flash4.0 (V3)		✓		
C13949-50U	Global Shutter CMOS Board Camera (12M)		✓		
C13770-50U	Global Shutter CMOS Board Camera (5M)		✓		
C13752-50U	Global Shutter CMOS Board Camera (3M)		✓		
C14041-10U	InGaAs QVGA Camera		✓		
C14041-20U	InGaAs QVGA Camera		✓		
C12741-03	InGaAs VGA Camera		✓		
C11440-62U	ORCA-Flash4.0 Board		✓		
C11440-52U	ORCA-Flash4.0 Board		✓		
C11440-52U30					
C11440-42U	ORCA-Flash4.0 LT		✓		
C11440-42U30					
C11440-36U	Global Shutter CMOS Camera		✓		
C11440-22CU	ORCA-Flash4.0 (V2)		✓		
C14300	X-Ray		✓		
C11800	X-Ray Line Dual Energy		✓		
C12849-111U	X-Ray CMOS		✓		
C12849-112U					
C12849-101U	X-Ray CMOS		✓		
C12849-102U					
C10400	X-Ray	✓		Windows 10 (*25) 32-bit / 64-bit (x64)	
C10650	X-Ray TDI	✓			
C10990	CCD Board camera	✓			
C10633-13/23	InGaAs camera	✓			
C9728DK-10	Flat panel sensor	✓		Windows 7 / 8.1 32-bit / 64-bit (x64)	
C9730DK-10		✓			
C9732DK-11		✓			

1394 OHCI

DCAM Versions

DCAM Module	19.12.322.5902	(for 32-bit)
	19.12.642.5902	(for 64-bit)
DRIVER	10.0.0.5902	

Cameras

Cameras	Nickname	400Mbps Max Card Speed	800Mbps Max Card Speed	Support OS	Note
C9100-24B	ImagEM X2 1K	n/a	BEST	Windows 10 (*25) 32-bit / 64-bit (x64)	
C11090-22B	ORCA-II	n/a	BEST		
C9100-23B	ImagEM X2	n/a	BEST	Windowsf 8.1 32-bit / 64-bit (x64)	
C10600-10B	ORCA-R2	poor	BEST		(*1)
C4742-80-12AG	ORCA-AG	good	BEST	Windows 7 (*20)/(*21) 32-bit / 64-bit (x64)	
C8484-xxG02	ORCA-xxG	good	BEST		
C9664-01G02		good	BEST		

Required

The IEEE-1394 card must have OHCI compatibility with Microsoft's Inbox Drivers.

Recommendations

- PCI Express x1 cards are better than PCI 32-bit/64-bit cards in desktops.
- ExpressCard is better than CardBus and PCMCIA cards in notebooks.
- Avago Technologies / LSI FW643 is the best PHY/Link IC for an IEEE-1394 interface card.
- It is highly recommended to disable C-state processor control in your PC's BIOS, else you may get sporadic corrupted images transferred to the PC. See Note (*23).

Notes

*New: New supported hardware or OS from the 19.10.5847 release.

*1: C10600-10B (ORCA-R2) and C11254-10B(ORCA-D2) have limitations for full performance with IEEE-1394 400Mbps port.

*2: This note is deprecated.

*3: This note is deprecated.

*4: This note is deprecated.

*5: This note is deprecated.

*6: This note is deprecated.

*7: This note is deprecated.

*8: This note is deprecated.

*9: This note is deprecated.

*10: This note is deprecated.

*11: This note is deprecated.

*12: This note is deprecated.

*13: This note is deprecated.

*14: This note is deprecated.

*15: PCIe x4 and x8 slot compatibility - For example, this is information comparing PCIe slot compatibility for some known Dell Precision Workstations and Dell Optiplex Desktops:

>Workstation T3600 or newer series

All slots are clearly labeled on the motherboard with their maximum link negotiation speed, width, as well as max power.

>Workstation T7500

Slot 2 PCI Express Gen2 x16 with support for up to 300W, full length, full height. Can support x8, x4 and x1 down shift.
(All Class Code cards should work, but certain cards may fall outside of spec. Report any issues.)

Slot 4 PCI Express Gen2 x16 with support for up to 225W, full length, full height. Can support x8, x4 and x1 down shift.
(All Class Code cards should work, but certain cards may fall outside of spec. Report any issues.)

Slot 1 PCI Express Gen2 x16 wired as x8 slot with support for up to 75W – half length, full height.
Can support x8, x4 and x1 down shift.

Slot 3 PCI Express Gen2 x16 wired as x8 slot with support for up to 75W – full length, full height.
Can support x8, x4 and x1 down shift.

Slot 6 PCI Express Gen2 x16 wired as x4 slot with support for up to 75W - full length, full height.
Can support x4, x1 down shift.

>Workstation T5500

Slots 2,4 PCI Express x16 with support for up to 225W – full length, full height. Can support x8, x4 and x1 down shift.
(All Class Code cards should work, but certain cards may fall outside of spec. Report any issues.)

Slot 1 PCI Express Gen2 x16 wired as x8 slot with support for up to 75W – half length, full height.
Can support x8, x4 and x1 down shift.

Slot 3 PCI Express Gen2 x16 wired as x8 slot with support for up to 75W – full length, full height.
Can support x8, x4 and x1 down shift.

>Workstation T3500

Slots 2,4 PCI Express x16 with support for up to 225W – full length, full height. Can support x8, x4 and x1 down shift.
(All Class Code cards should work, but certain cards may fall outside of spec. Report any issues.)

Slots 1,3 PCI Express x8 wired as x4 slot with support for up to 75W – half length, full height.
Can support x4 and x1 down shift.

>Workstation T1500

Slot 1 PCI Express x16 with support for up to 75W – half length, full height. Can support x8, x4, and x1 down shift.
(All Class Code cards should work, but certain cards may fall outside of spec. Report any issues.)

Here is some compatibility information for some known HP Workstations:

> HP Z220, Z420, Z620, and Z820 Workstations

- View the [Maintenance and Service Guide](#)
 - Component replacement guidelines
 - Expansion slots

> HP Z210 CMT Workstation

- View the [Maintenance and Service Guide](#)
 - Component replacement guidelines
 - Expansion slots

Notes – Cont'd

>Z400 Workstation (Taken from the HP Z400 Maintenance and Service Guide)


Expansion card slot identification

This section identifies and describes workstation expansion card slots, and presents card configuration information.

Slot identification and description

The following figure identifies workstation expansion card slots.

Identifying expansion card slots

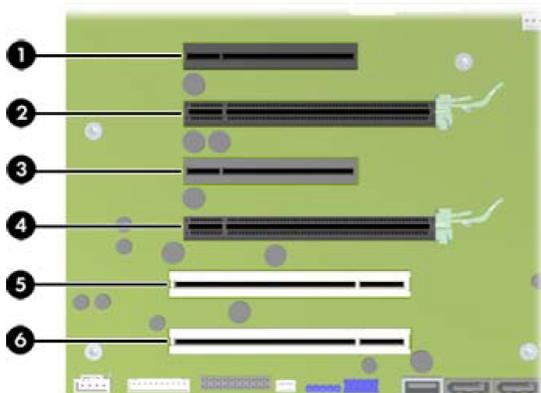

The following table describes the workstation PCIe card slots.

Slot	Type	Slot power (Maximum)
1	PCIe2 - x8(4)	25W
2	PCIe2 - x16	75W
3	PCIe - x8(4)	25W
4	PCIe2 - x16	75W

NOTE: The x1, x4, x8, and x16 designators describe the mechanical length of the slot. The number in parentheses shows how many electrical PCIe lanes are routed to the expansion slot. For example, x16(8) means that the expansion slot is mechanically a x16 length connector, with eight PCIe lanes connected. A x16 PCIe card runs at the bandwidth of the slot it is plugged into.

Slots one and three use open-ended PCIe connectors, so a PCIe x16 card can be inserted. Graphics cards greater than 75 watts require the use of an auxiliary power cable adapter.

Expansion card

PCIe I/O slots can support other PCIe cards with less bus bandwidth than what is physically defined for the slot.

Expansion card slot description

NOTE: The x1, x4, x8, and x16 designators describe the mechanical length of the slot. The number in parentheses shows how many electrical PCIe lanes are routed to the expansion slot. For example, x16(8) means that the expansion slot is mechanically a x16 length connector, with eight PCIe lanes connected. A x16 PCIe card runs at the bandwidth of the slot it is plugged into.

Use the following table to determine PCIe card compatibility.

Workstation PCIe compatibility matrix

Slot	Mechanical compatibility	Electrical compatibility
1	x8 connector, open-ended*	PCIe2 (x1, x4)
2	x16 connector	PCIe2 (x1, x4, x8, x16)
3	x8 connector, open-ended*	PCIe (x1, x4)
4	x16 connector	PCIe2 (x1, x4, x8, x16)

*This slot is open-ended so a PCIe x16 card can be inserted.


- Slots one, two, and four are PCIe GEN2 slots.
- Graphics cards greater than 75 watts require the use of an auxiliary power cable adapter.

>Z800 Workstation (Taken from the HP Z800 Maintenance and Service Guide)

Expansion card slot description

The following figure identifies workstation expansion card slots.

Identifying expansion card slots

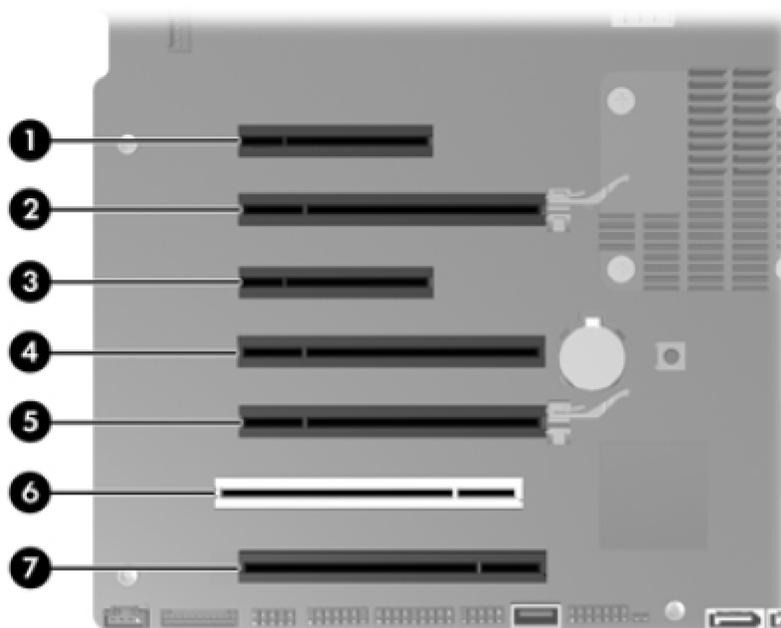

The following table describes the workstation expansion card slots.

Expansion card slot description and compatibility

Slot	Type	Mechanical compatibility	Electrical compatibility	Slot power (Maximum)
1 ¹	PCIe2 x8(4) open-ended	x1, x4, x8, x16	x1, x4	25W
2 ²	PCIe2 x16 ³	x1, x4, x8, x16	x1, x4, x8, x16	75W
3	PCIe x8(4) open-ended	x1, x4, x8, x16	x1, x4	25W
4 ¹	PCIe2 x16(8)	x1, x4, x8, x16	x1, x4, x8	25W
5 ¹	PCIe2 x16 ⁴	x1, x4, x8, x16	x1, x4, x8, x16	75W
7	PCIe2 x16(8)	x1, x4, x8, x16	x1, x4, x8	25W

¹ PCIe GEN2 slot.

² Primary graphics slot

⁴ Secondary graphics slot

NOTE: The x1, x4, x8, and x16 designators describe the number of electrical PCIe lanes routed to an expansion slot. For example, x16(8) means that the expansion slot is mechanically a x16 length connector, with eight PCIe lanes connected.


Slots one and three use open-ended PCIe connectors, so a PCIe x16 card can be inserted. This allows the workstation to support more PCIe x16 expansion cards. A x16 card typically trains and runs at the maximum lane width available by the expansion slot. The card runs at the reduced lane width, with a corresponding performance decrease.

A x16 graphics card runs at the bandwidth of the slot it is plugged into. The slot bandwidth can be x1, x4, x8, or x16. A PCIe card can be plugged into a slot with a lesser speed. It runs at that speed with a commensurate performance decrease. (Older graphics cards may not comply with this.)

- *16: This note is deprecated.
- *17: This note is deprecated.
- *18: The C11440-10C (ORCA-Flash 2.8) with a Phoenix PE1 card requires the camera firmware version to be 1.1 or newer.
- *19: These hardware must be connected to a USB 3.0 compliant bus. If these hardware are connected to a [Renesas \$\mu\$ PD720202](https://www.renesas.com/en/products/usb-controllers/usb-3.0-host-controllers/usb-3.0-host-controller-drivers/30230) chipset controller in **Windows 7**, download this Renesas driver:

ftp://60.248.38.84/cat_106/30230_dr.zip

- (a) After you download this driver set, extract the contents to a folder on your PC.
- (b) Go to Add or remove programs, scroll down and look for this entry, and if you see it, Uninstall it:

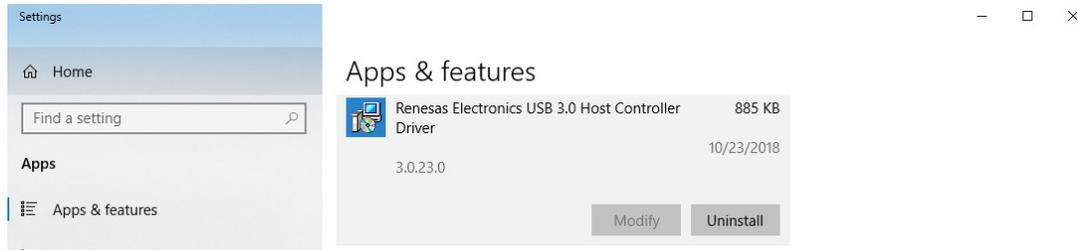

- (c) Install this driver set with this exact command line (including switches) from within the folder you extracted the driver set:

```
RENESAS-30230-setup.exe /s /v"NO_MONITOR=1 NO_UTILITY=1 SET_SELSUS_MODE1=1
U1U2_DISABLE=1 U1U2_UTL_DISABLE=1 U1U2_PROPERTY_DISABLE=1"
```

This exact command line can be downloaded from this link:

<ftp://hamamatsu.hostedftp.com/~hamamatsu/DCAM-API/Renesas/RENESAS-30230-setup.cmd>

If you download this .CMD file to the same folder as RENESAS-30230-setup.exe from step (a), you should be able to execute it AsAdmin to run with the exact noted command line with switches to install the driver properly.

- *20 For 1394 and Windows 7 only, it is recommended that you install Service Pack 1:

<http://support.microsoft.com/kb/976932>

Microsoft has made many improvements to their 1394 driver stack which are incorporated in SP1 and not necessarily documented or included in RTM hotfixes.

- *21 For 1394 and Windows 7 only, if your experiment requires you to make many captures with various binning, sub-array, scan speed, and/or data type settings over hours of operation (usually more than 1 hour), there is a known bug inside Microsoft's latest driver standard driver stack which could present itself through our API to the host application as an "Out of Resource" or "No Resource" error. When this happens, there is no way to recover capturing again from our 1394 device unless you exit your current application, power cycle the device, and restart the experiment manually. We have traced the root cause of the problem. It happens when our lower drivers cannot free 1394 Isochronous bandwidth properly through the Microsoft drivers, we can no longer re-use that bandwidth, and further allocations we require will fail.

If you experience the above known issue, it is recommended you install SP1 if you are not at this level:

<http://support.microsoft.com/kb/976932>


Microsoft has released an official hotfix to address this issue:

<http://support.microsoft.com/kb/2524249>


If you configured your IEEE1394 Bus controller driver for Legacy per a previous Compatibility Note recommendation, you should revert your IEEE 1394 Bus Controller driver away from Legacy and back to the Standard driver to use this hotfix and be the most stable. The Legacy driver is provided in Windows 7 only for backward driver compatibility. It will likely not evolve in stability and function beyond its current state.

- *22 The FireBird 2PE8 cards operate optimally with PCIe x8 Gen2 slots. Please make sure you install this card into an electrically compatible PCIe x8 Gen2 slot. If you have a Dell or HP Workstation, please reference to (*15) above.
- *23 For all PCIe interface options, it is highly recommended to disable C-state processor control in your PC's BIOS. Most PC manufacturers (such as Dell and HP) ship with C-state processor control enabled from the factory. When this is enabled, the processor may drop in and out of S0 maximum power state – when this happens, if you are transferring image data across the PCIe bus, it may get interrupted to the point where data becomes corrupted or lost. It is also known that C-state being enabled can affect the overall maximum bandwidth of particular PCIe slots on the motherboard. For critical image capturing, insure S0 power state is always maximum – the best way to do it is to tweak the BIOS for such control.

For example, if you have a Dell T3500, T5500, and/or T7500, the setting appears in the BIOS under Performance:

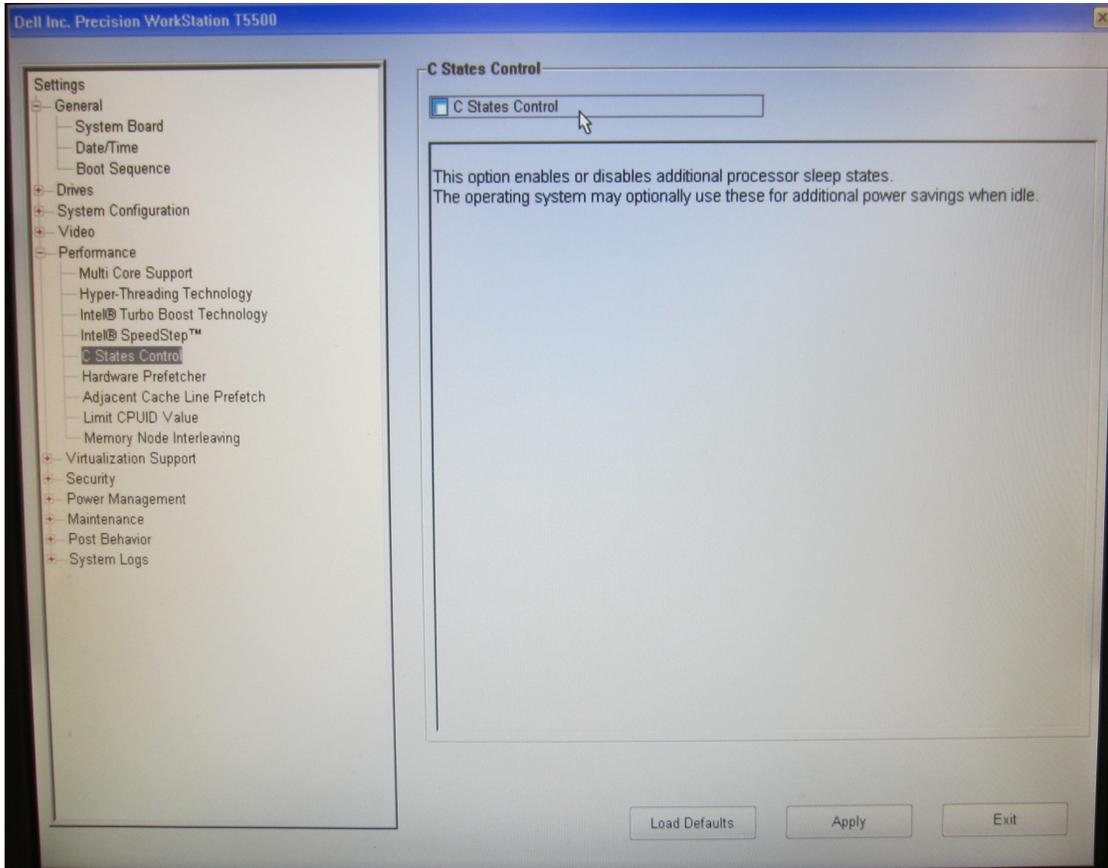

If you have an HP Z400 or Z800 Workstation, it appears in the BIOS here:

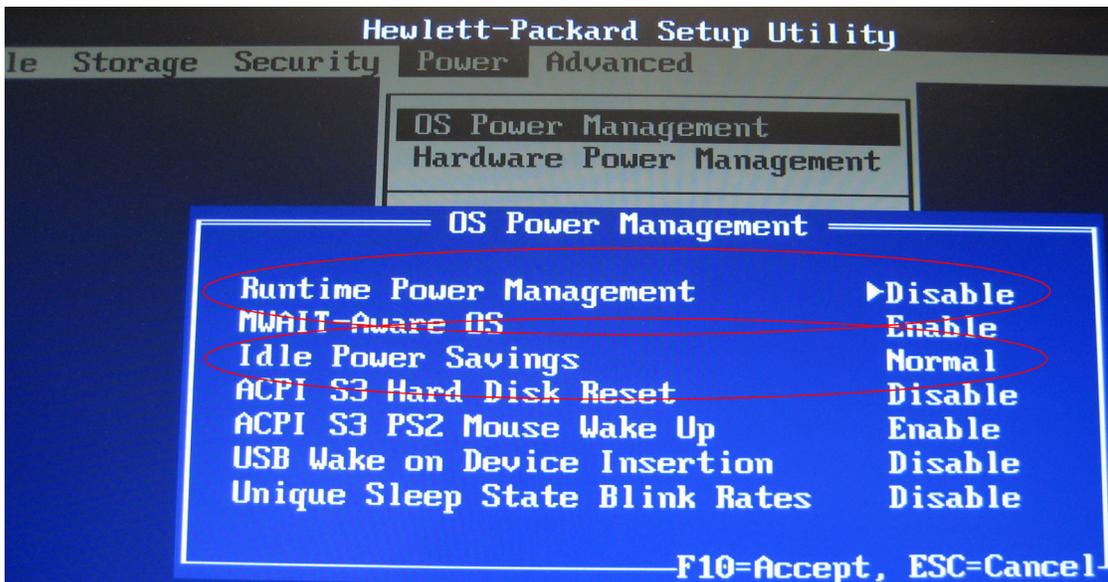


Set Runtime Power Management to Disable and Idle Power Savings to Normal. For a further description from HP (Taken from Computer Setup(F10) Utility):

Computer Setup—Power	
Option	Description
OS Power Management	<ul style="list-style-type: none"> Runtime Power Management— Enable/Disable. Allows certain operating systems to reduce processor voltage and frequency when the current software load does not require the full capabilities of the processor. Idle Power Savings—Extended/Normal. Allows certain operating systems to decrease the processors power consumption when the processor is idle.

- *24 These hardware must be connected to a USB 3.0 compliant bus, and the drivers for the USB 3.0 chipset controller must be operational in Device Manager (check the card or PC manufacturer's website for proper drivers if Windows 7. Windows 8 and higher ship with in-box USB xHCI compliant drivers from Microsoft that can work with any known USB3 chipset controller), and the USB 3.0 chipset controller must be installed/operating on a PCI Express / ExpressCard **Gen2** (5GT/s) capable slot / bus. [Renesas \$\mu\$ PD720202](#) is the most compatible USB host controller chipset with these hardware. When using the Renesas μ PD720202 controller and these hardware, download this Renesas driver, **even if you are using Windows 8 or higher** – the Microsoft in-box drivers do not work as well as the Renesas drivers when coupled with these hardware:

ftp://60.248.38.84/cat_106/30230_dr.zip

- (a) After you download this driver set, extract the contents to a folder on your PC.
 (b) Go to Add or remove programs, scroll down and look for this entry, and if you see it, Uninstall it:

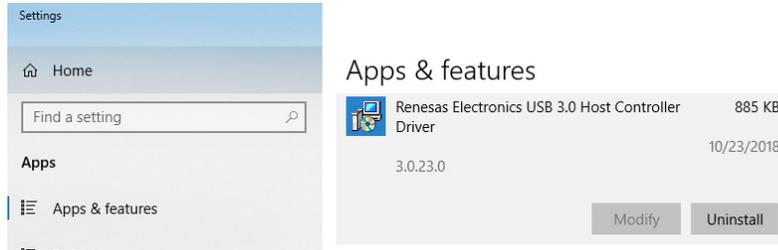

- (c) Install this driver set with this exact command line (including switches) from within the folder you extracted the driver set:


```
RENESAS-30230-setup.exe /s /v"NO_MONITOR=1 NO_UTILITY=1 SET_SELSUS_MODE1=1
U1U2_DISABLE=1 U1U2_UTL_DISABLE=1 U1U2_PROPERTY_DISABLE=1"
```

This exact command line can be downloaded from this link:

<ftp://hamamatsu.hostedftp.com/~hamamatsu/DCAM-API/Renesas/RENESAS-30230-setup.cmd>

If you download this .CMD file to the same folder as RENESAS-30230-setup.exe from step (a), you should be able to execute it AsAdmin to run with the exact noted command line with switches to install the driver properly.

If you install the Renesas driver to Windows 8 or higher, it will change the driver for every Renesas μ PD720201/2 controller present in your system during the time of installation. If you have multiple USB 3.0 devices (ex – a C13440-20CU ORCA-Flash4.0 (V3)) in the same system connected to independent Renesas μ PD720202 controllers than the hardware mentioned above, that hardware may work with better performance if the Microsoft in-box drivers are used. For those other USB 3.0 devices, you can use Device Manager with View devices by connection, then Update Driver to the Renesas controller connected to the other devices manually to switch those controllers back to the Microsoft in-box driver. Device Manager could then look like this example with the Renesas driver for one controller, and the Microsoft in-box driver for the other:

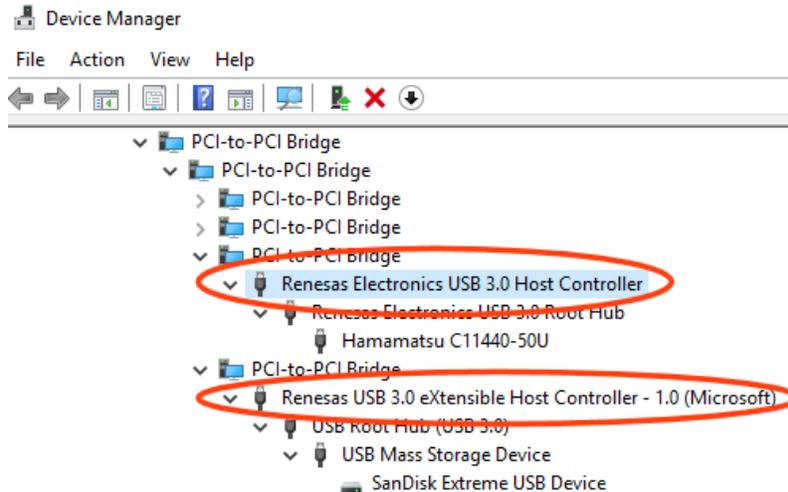

*25 Verified using Windows10 Ver1903. If you are using Windows 10 OS newer than Ver1903, it may not work correctly.