

Mini-spectrometer

[TF series]

C14486GA



For near-infrared range, compact/thin case with InGaAs linear image sensor mounted

The mini-spectrometer TF series is a polychromator provided in a compact, thin case that houses optical elements, image sensor, and driver circuit. Spectrum data can be acquired by guiding measurement light into a mini-spectrometer through an optical fiber and transferring the measured results to a PC via the USB connection. Moreover, the trigger function that can be also used for short-term integration enables spectroscopic measurement of pulse emissions.

The product includes free evaluation software with functions for setting measurement conditions, acquiring and saving data, drawing graphs, and so on. Furthermore, the DLL function specifications are disclosed, so users can create their original measurement software programs.

Features

- Compact, thin case
- InGaAs linear image sensor built-in
- With trigger function
- High throughput using quartz transmission grating
- External power supply is not necessary (USB bus powered).
- Easy installation in devices
- Stores wavelength conversion factor*1 in internal memory

*1: A conversion factor for converting the image sensor pixel number into a wavelength. A calculation factor for converting the A/D converted count into the input light level is not provided.

Applications

- Sugar content of foods, moisture measurement
- Film thickness measurement

Structure

Parameter	Specification	Units
Dimensions (W × D × H)	80 × 60 × 12	mm
Weight	88	g
Image sensor	InGaAs linear image sensor	-
Number of pixels	256	pixels
Slit*2 (H × V)	25 × 250	μm
NA*3	0.22	-
Connector for optical fiber	SMA905D	-

*2: Input slit aperture size

*3: Numeric aperture (solid angle)

Absolute maximum ratings

Parameter	Value	Units
Operating temperature*4	+5 to +50	°C
Storage temperature*4	-20 to +70	°C

*4: No dew condensation

When there is a temperature difference between a product and the surrounding area in high humidity environment, dew condensation may occur on the product surface. Dew condensation on the product may cause deterioration in characteristics and reliability.

Note: Exceeding the absolute maximum ratings even momentarily may cause a drop in product quality. Always be sure to use the product within the absolute maximum ratings.

Optical characteristics

Parameter	Specification	Units
Spectral response range	950 to 1700	nm
Spectral resolution (half width) ^{*5}	Typ.	5.0
	Max.	7.0
Wavelength reproducibility ^{*6}	-0.4 to +0.4	nm
Wavelength temperature dependence	-0.05 to +0.05	nm/°C
Spectral stray light ^{*5 *7}	-33 max.	dB

*5: When the slit in the table in "Structure" is used. The spectral resolution depends on the slit.


*6: Measured under constant light input conditions

*7: The ratio of the count measured when a 1300 nm light is input to the count measured when a 1300 ± 40 nm light is input.

Electrical characteristics


Parameter	Specification	Unit
A/D conversion	16	bit
Integration time	1 to 100000	μs
Interface	USB 2.0	-
USB bus power current consumption	Typ.	220
	Max.	250

Spectral response (typical example)

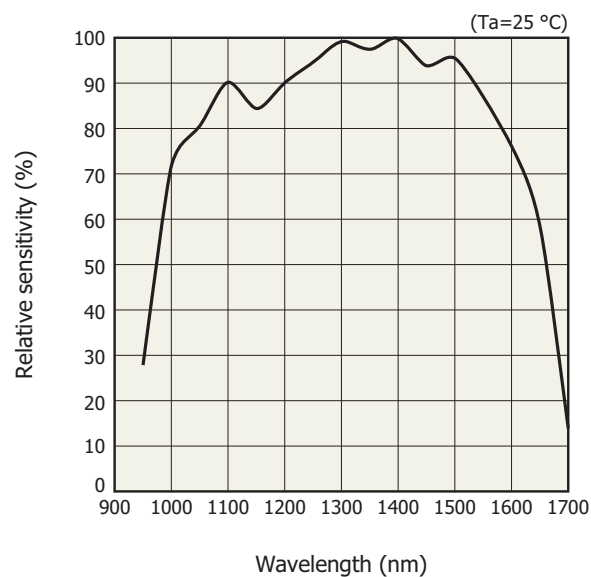


KACCB0509EA

Spectral resolution vs. wavelength (typical example)

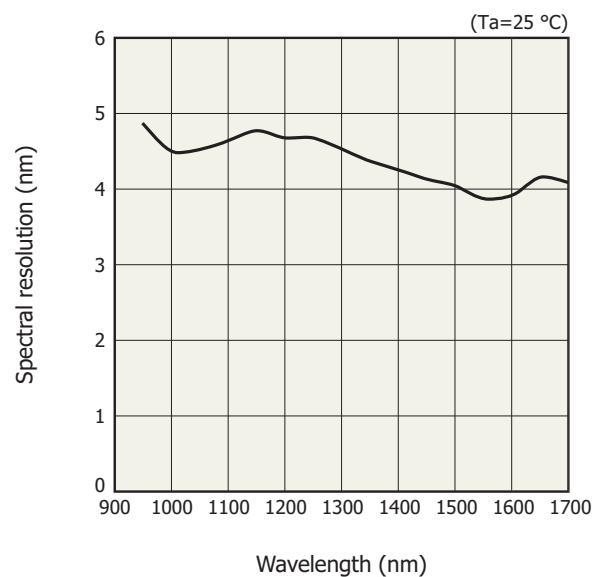

KACCB0510EA


Linearity (typical example)

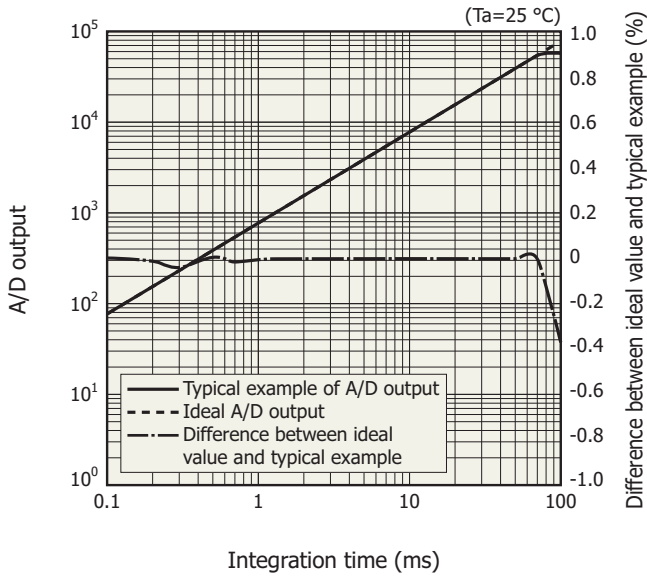

KACCB0511EA

A/D output is the output with dark output subtracted when light is input. The difference between the ideal value and typical example contains a measurement error. The smaller the A/D output, the larger the measurement error

Dark output vs. temperature (typical example)

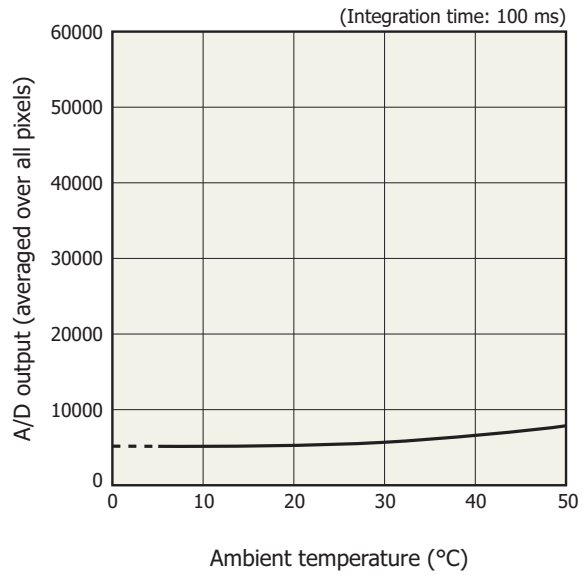


KACCB0512EA

A/D output is the sum of the sensor and circuit offset outputs and the sensor dark output.

Optical component layout


The mini-spectrometer TF series employs a transmission holographic grating made of quartz and an optical system arranged on a robust optical base to produce high throughput and highly accurate optical characteristics.

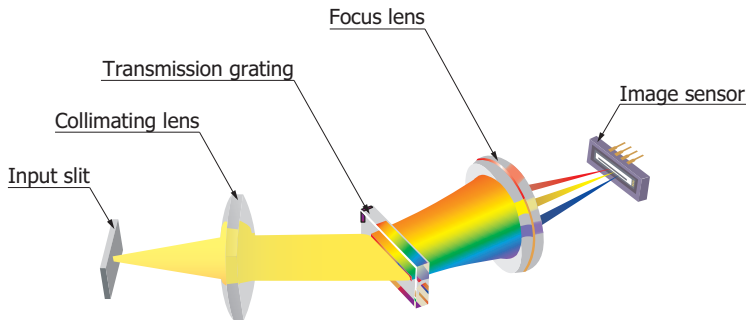

KACCC0917EA

Connection example (transmitted light measurement)

Spectrum data can be acquired by guiding measurement light into a mini-spectrometer through an optical fiber and transferring the measured results to a PC via the USB connection. Since there are no moving parts inside the device, constantly stable measurements can be expected. Moreover, the optical guiding section uses an optical fiber making connection to the measured object flexible.

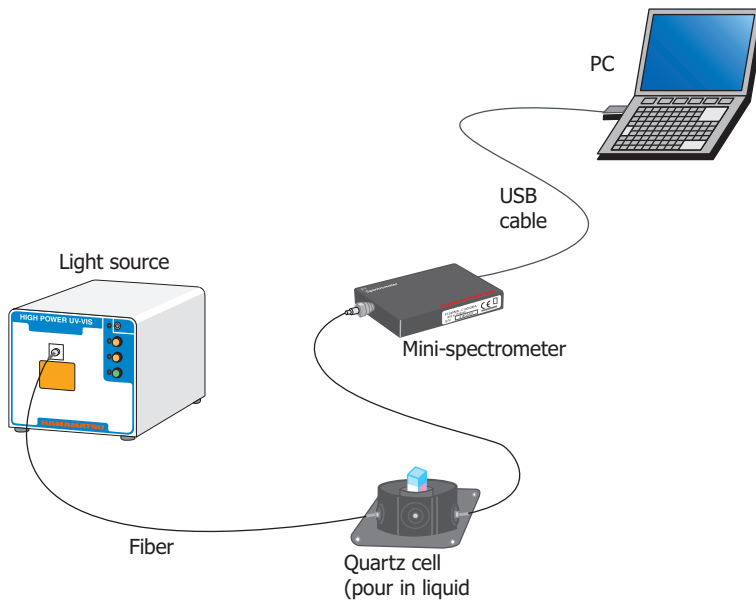


KACCC0767EB

Trigger operation modes

In the C14486GA, the following trigger operation modes are available. You can switch between these modes from the evaluation software supplied with the C14486GA.

(1) Asynchronous data measurement at software trigger input


The first piece of digital data that is converted after a software trigger is applied from the PC is acquired.

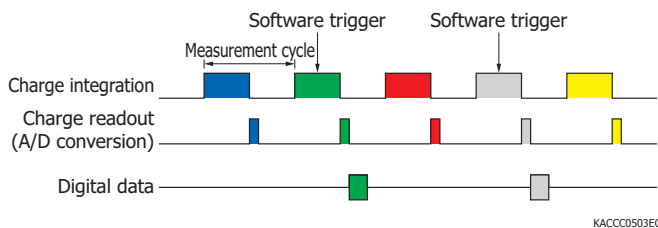

KACCC0503EC

(2) Synchronous data measurement at software trigger input


Sensor operation (integration) starts when a software trigger is applied from the PC.

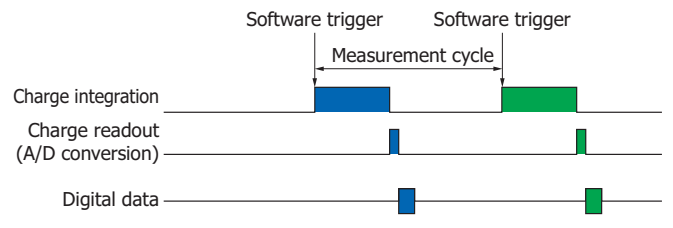

KACCC0505EB


(3) Asynchronous data measurement at external trigger input

The first piece of digital data that is converted after an external trigger edge (rising or falling edge can be specified) is applied to the external trigger terminal is acquired.

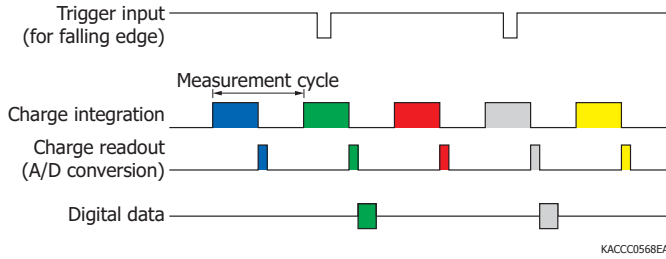


(4) Synchronous data measurement at external trigger input

Sensor operation (integration) starts when an external trigger edge (rising or falling edge can be specified) is applied to the external trigger terminal, and then the digital data is acquired.

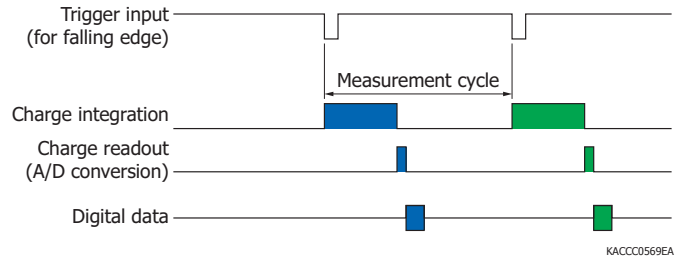


(5) Asynchronous data measurement at external trigger input level

Digital data is acquired when an external trigger (high level or low level can be specified) is applied to the external trigger terminal.

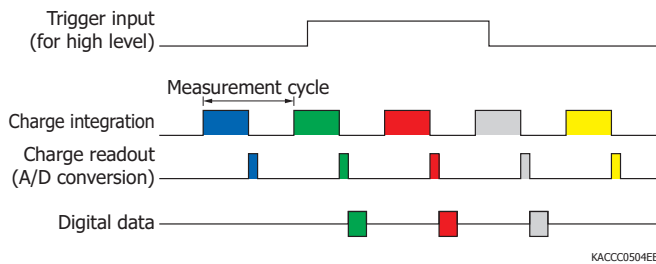

(6) Synchronous data measurement at external trigger input level


Sensor operation (integration) starts when a trigger (high level or low level can be specified) is applied to the external trigger terminal, and then the digital data is acquired.

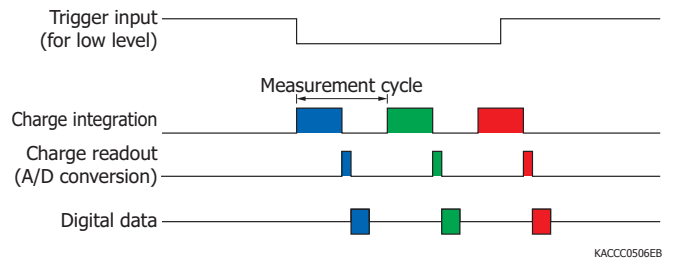

In any of the modes (1) to (6), if the trigger input cycle is shorter than the measurement cycle of the spectrometer, the input trigger is ignored.

(7) External trigger signal output


The start timing (pulse width: 10 μ s) of integration can be output from the external trigger terminal (trigger output edge: rising or falling edge can be specified).

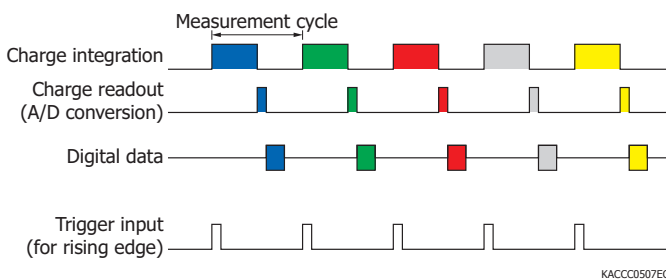

❏ Evaluation software (accessory)

By installing the evaluation software (SpecEvaluationUSB2.exe)*8 into a PC, you can perform the following basic operations.

- Acquire and save measured data
- Set measurement conditions
- Module information acquisition (wavelength conversion factor, mini-spectrometer type, etc.)
- Display graphs
- Arithmetic functions
 - Pixel number to wavelength conversion
 - Calculation in comparison with reference data (transmittance, reflectance)
 - Dark subtraction
 - Gaussian approximation (peak position and count, FWHM)

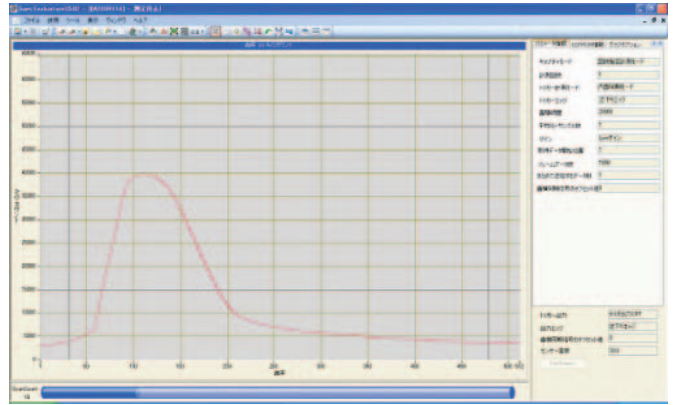

Note: Up to eight mini-spectrometers can be connected to a single PC and used.

*8: Compatible OS

- Microsoft® Windows® 7 (32-bit, 64-bit)
- Microsoft Windows 8 (32-bit, 64-bit)
- Microsoft Windows 10 Professional (32-bit, 64-bit)


A DLL for controlling the hardware is available.

Users can develop original measurement programs using the following development platform.

- Microsoft Visual Studio® 2008 (SP1) Visual C++®
- Microsoft Visual Studio 2008 (SP1) Visual Basic®

Note: Microsoft, Windows, Visual Studio, Visual C++, and Visual Basic are either registered trademarks or trademarks of Microsoft Corporation in the United States and/or other countries.

❏ Dimensional outline (unit: mm)

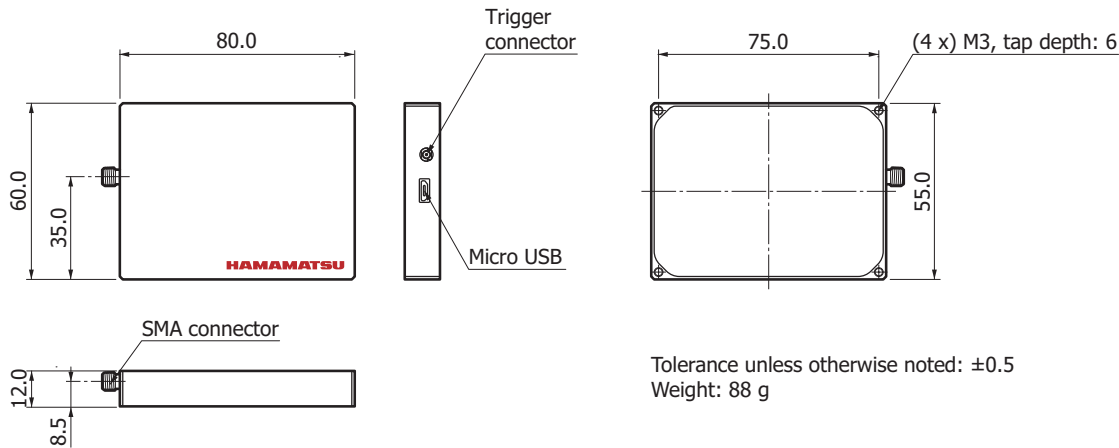

KACCA0415EA

Accessories

- USB cable
- Dedicated software (evaluation software, sample software, DLL)


Options (sold separately)

- Input optical fiber

Type no.	Product name	Core diameter (μm)	Specifications
A9763-01	Fiber for visible/near infrared range	600	NA=0.22, length=1.5 m, low cost With SMA905D connector on each end
A9763-05		400	NA=0.22, length=1.5 m, small bending radius at fiber section With SMA905D connector on each end

- External trigger input coaxial cable A12763

Dimensional outline (unit: mm)

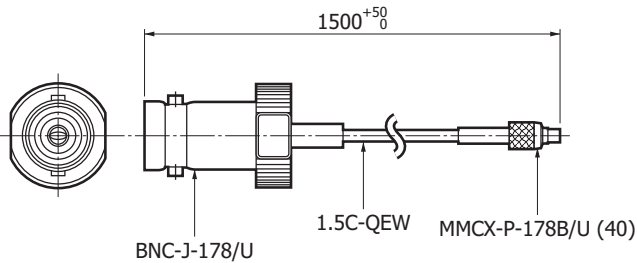

Mini-spectrometer lineup

Type no.	Type	Spectral response range(nm)													Spectral resolution max. (nm)	Image sensor		
		200	400	600	800	1000	1200	1400	1600	1800	2000	2200	2400	2600				
C10082CA	TM-UV/VIS-CCD High sensitivity																6	Back-illuminated type CCD Image sensor
C10082CAH	TM-UV/VIS-CCD High resolution		200 to 800														1*	
C10082MD	TM-UV/VIS-MOS Wide dynamic range																6	CMOS linear image sensor
C10083CA	TM-VIS/NIR-CCD High sensitivity																8 (Wavelength 320 to 900 nm)	Back-thinned CCD image sensor
C10083CAH	TM-VIS/NIR-CCD High resolution																1* (Wavelength 320 to 900 nm)	
C10083MD	TM-VIS/NIR-MOS Wide dynamic range																8	CMOS linear image sensor
C11697MB	TM-VIS/NIR-MOS-II Trigger compatibility																8	High-sensitivity CMOS linear image sensor
C9404CA	TG-UV-CCD High sensitivity		200 to 400														3	Back-thinned CCD image sensor
C9404CAH	TG-UV-CCD High resolution																1*	Back-thinned CCD image sensor
C9405CB	TG-SWNIR-CCD-II IR enhanced																5 (Wavelength 550 to 900 nm)	IR-enhanced back-thinned CCD image sensor
C11713CA	TG-RAMAN-I High resolution																0.3*	Back-thinned CCD image sensor
C11714CB	TG-RAMAN-II High resolution																0.3*	IR-enhanced back-thinned CCD image sensor
C11482GA	TG2-NIR Non-cooled type																7	InGaAs linear image sensor
C9913GC	TG-cooled NIR-I Low noise(Cooled type)																7	
C9914GB	TG-cooled NIR-II Low noise(Cooled type)																8	
C11118GA	TG-cooled NIR-III Low noise(Cooled type)																20	
C13555MA	TF-VIS-MOS-II Compact, thin case																3	High-sensitivity CMOS linear image sensor
C13053MA	TF-SWIR-MOS-II Compact, thin case																3.5	
C14486GA	TF-NIR Compact, thin case																5*	InGaAs linear image sensor
C13054MA	TF-RAMAN Compact, thin case																0.4*	High-sensitivity CMOS linear image sensor
C14214MA	TF-RAMAN Compact, thin case																0.6	
C11007MA	RC-VIS-MOS Spectrometer module																9	CMOS linear image sensor
C11008MA	RC-SWNIR-MOS Spectrometer module																8	IR-enhanced CMOS linear image sensor

* Typ. Value

For incorporation into mobile measurement devices

Type no.	Type	Spectral response range(nm)													Spectral resolution max. (nm)	Image sensor		
		200	400	600	800	1000	1200	1400	1600	1800	2000	2200	2400	2600				
C11009MA	RC-VIS-MOS Spectrometer head																9	CMOS linear image sensor
C11010MA	RC-SWNIR-MOS Spectrometer head																8	IR-enhanced CMOS linear image sensor

For incorporation into mobile measurement devices (Ultra-compact)

Type no.	Type	Spectral response range(nm)													Spectral resolution max. (nm)	Image sensor		
		200	400	600	800	1000	1200	1400	1600	1800	2000	2200	2400	2600				
C11708MA	MS-SWNIR-MOS Spectrometer head																20	CMOS linear image sensor
C12666MA	Spectrometer head																15	CMOS linear image sensor
C12880MA	Spectrometer head																15	High-sensitivity CMOS linear image sensor

Related information

www.hamamatsu.com/sp/ssd/doc_en.html

■ Precautions

- Disclaimer
- Mini-spectrometers

■ Technical information

- Mini-spectrometers

Information described in this material is current as of August 2018.

Product specifications are subject to change without prior notice due to improvements or other reasons. This document has been carefully prepared and the information contained is believed to be accurate. In rare cases, however, there may be inaccuracies such as text errors. Before using these products, always contact us for the delivery specification sheet to check the latest specifications.

The product warranty is valid for one year after delivery and is limited to product repair or replacement for defects discovered and reported to us within that one year period. However, even if within the warranty period we accept absolutely no liability for any loss caused by natural disasters or improper product use. Copying or reprinting the contents described in this material in whole or in part is prohibited without our prior permission.

HAMAMATSU

www.hamamatsu.com

HAMAMATSU PHOTONICS K.K., Solid State Division

1126-1 Ichino-cho, Higashi-ku, Hamamatsu City, 435-8558 Japan, Telephone: (81) 53-434-3311, Fax: (81) 53-434-5184

U.S.A.: Hamamatsu Corporation: 360 Foothill Road, Bridgewater, N.J. 08807, U.S.A., Telephone: (1) 908-231-0960, Fax: (1) 908-231-1218, E-mail: usa@hamamatsu.com

Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49) 8152-375-0, Fax: (49) 8152-265-8, E-mail: info@hamamatsu.de

France: Hamamatsu Photonics France S.A.R.L.: 19, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: 33-(1) 69 53 71 00, Fax: 33-(1) 69 53 71 10, E-mail: infos@hamamatsu.fr

United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10 Tewin Road, Welwyn Garden City, Hertfordshire AL7 1BW, United Kingdom, Telephone: (44) 1707-294888, Fax: (44) 1707-325777, E-mail: info@hamamatsu.co.uk

North Europe: Hamamatsu Photonics Norden AB: Torshamnsgatan 35 16440 Kista, Sweden, Telephone: (46)8-509 031 00, Fax: (46)8-509 031 01, E-mail: info@hamamatsu.se

Italy: Hamamatsu Photonics Italia S.r.l.: Strada della Moia, 1 int. 6, 20020 Arese (Milano), Italy, Telephone: (39)02-93 58 17 33, Fax: (39)02-93 58 17 41, E-mail: info@hamamatsu.it

China: Hamamatsu Photonics (China) Co., Ltd.: B1201, Jiaming Center, No.27 Dongsanhuan Beilu, Chaoyang District, Beijing 100020, China, Telephone: (86) 10-6586-6006, Fax: (86) 10-6586-2866, E-mail: hpc@hamamatsu.com.cn

Taiwan: Hamamatsu Photonics Taiwan Co., Ltd.: 8F-3, No. 158, Section2, Gongdao 5th Road, East District, Hsinchu, 300, Taiwan R.O.C. Telephone: (886)03-659-0080, Fax: (886)03-659-0081, E-mail: info@hamamatsu.com.tw